





Introducing Proseek® Multiplex
Truly Scalable Immunoassays for
Biomarker Research

Protein Biomarkers

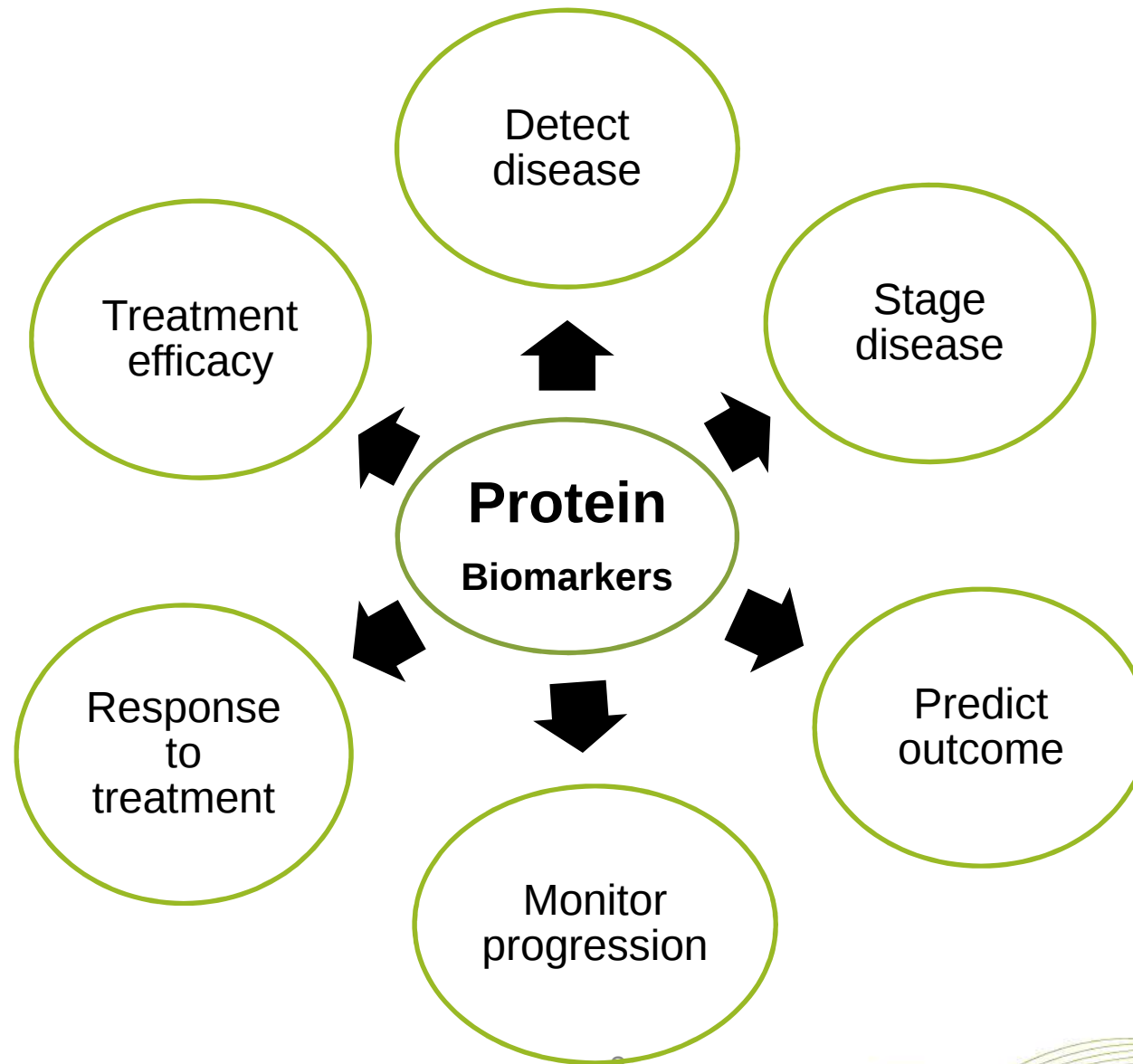


Development of new biomarkers

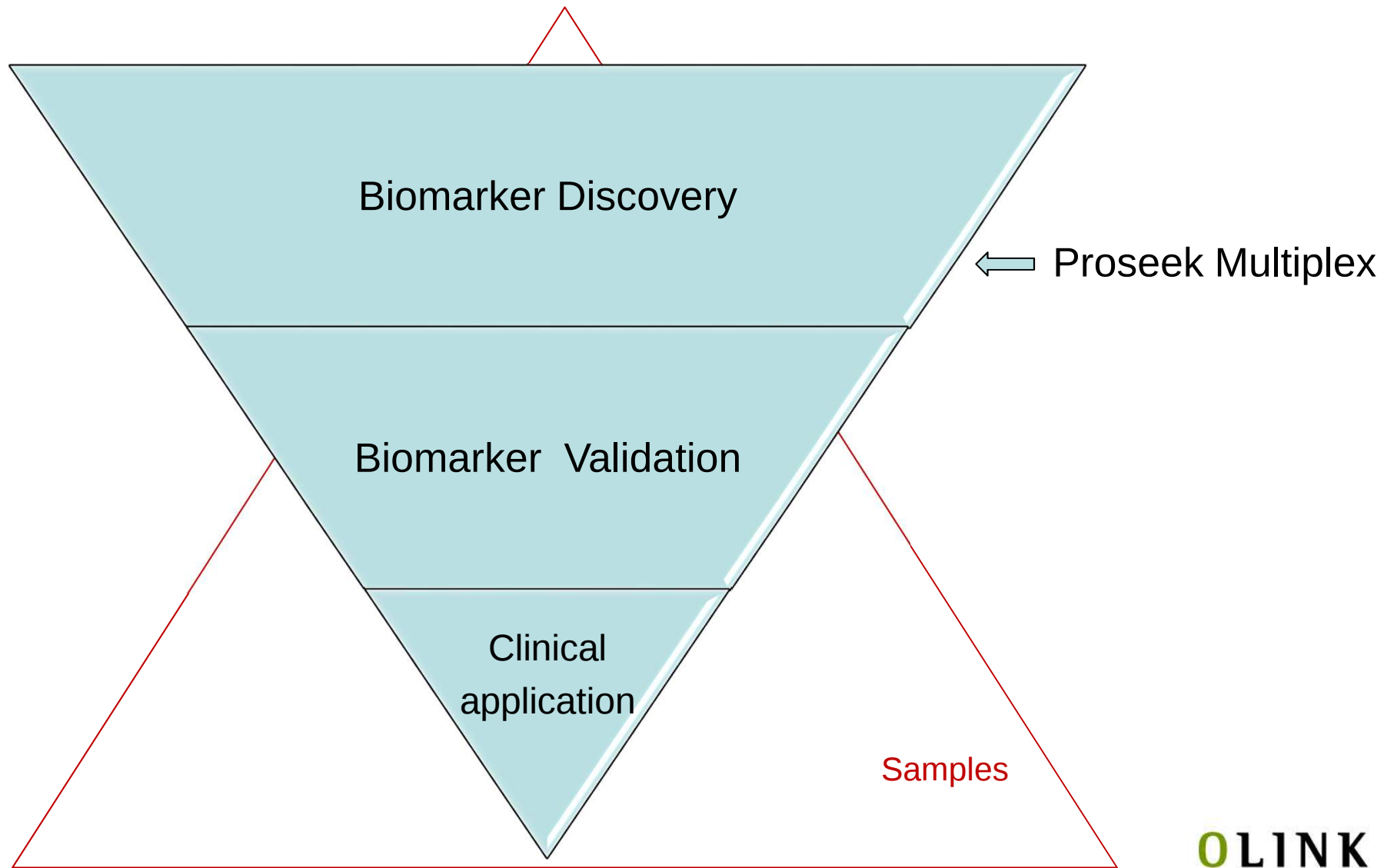

We address two of the major challenges

- ▶ High multiplexing: many analytes per sample : 92
- ▶ Low sample consumption : 1 uL

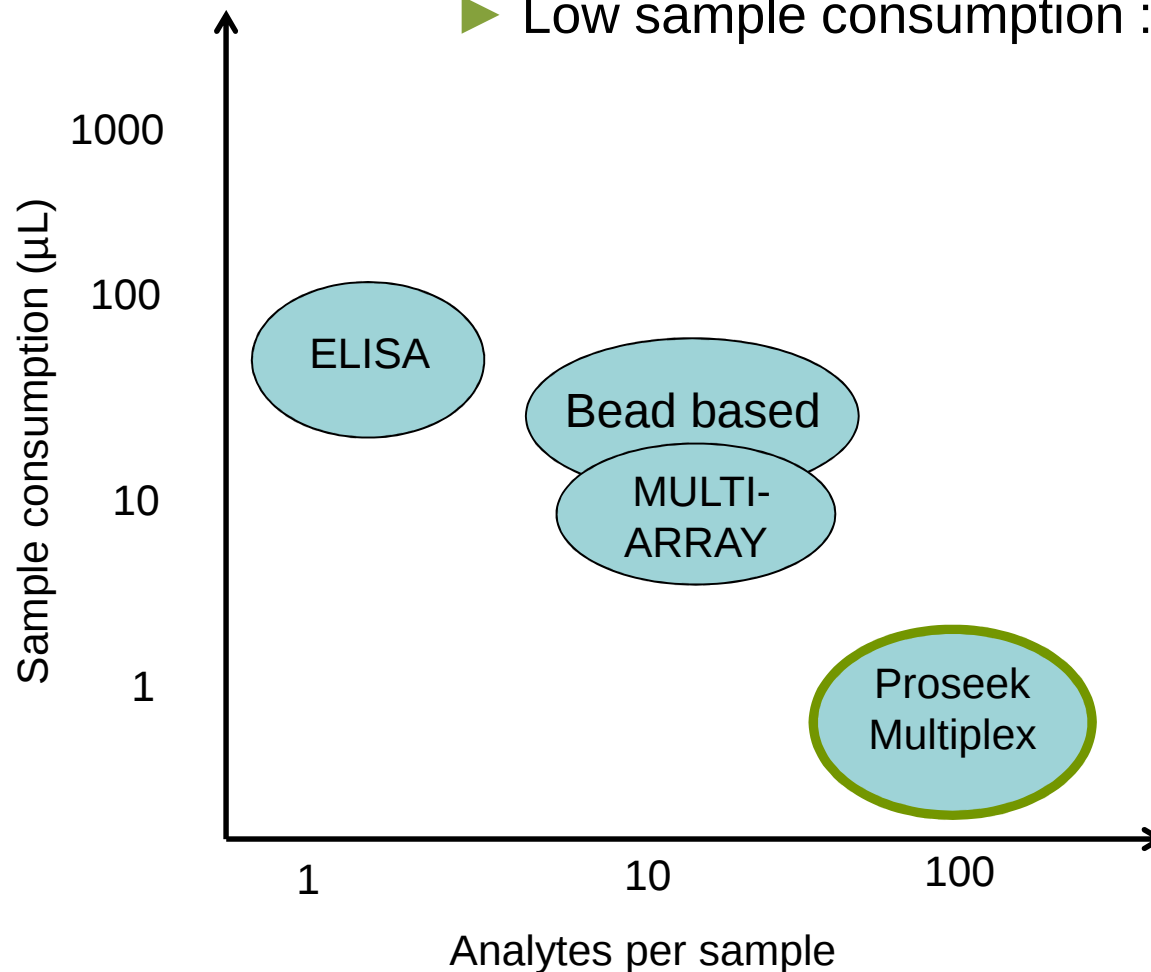

Why study many proteins at the same time?

- ▶ When you don't know exactly what you're looking for
 - Biomarker discovery
- ▶ Getting the complete picture
 - Proteins rarely act alone, for example in inflammation
- ▶ Discovering novel biomarker patterns for in vitro diagnostic use

Why study many proteins in small sample volumes?


- ▶ Enables access to precious biobanked serum and plasma
 - Retrospective biomarker research
 - From clinical trials, translational research
- ▶ Analysis of small tissue samples
 - Biopsies, even from needles
- ▶ Laboratory animals such as mice
 - Pharmacodynamics and toxicology studies

Why study many proteins in many samples?


- ▶ Getting statistical power in multivariate biomarker discovery requires many samples
- ▶ Correlating protein biomarker data to....
 - Multiple disease outcomes
 - Multiple therapeutic avenues
 - Multiple genetic blue prints (we are all different as individuals)

What does Proseek Multiplex do?

- ▶ Target multiple markers
 - 92 human proteins
- ▶ Small sample consumption
 - just one 1µL is required
- ▶ And most importantly, highly specific
 - Ensures the signal measured is from the correct protein !
 - Can not be done with competing technologies

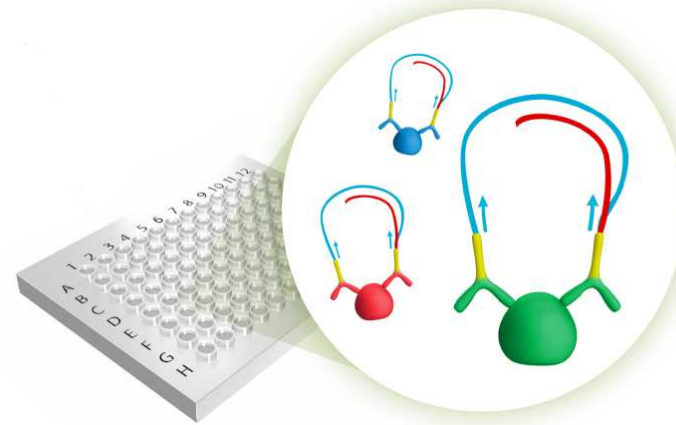

Countinous development of biomarker panels.....

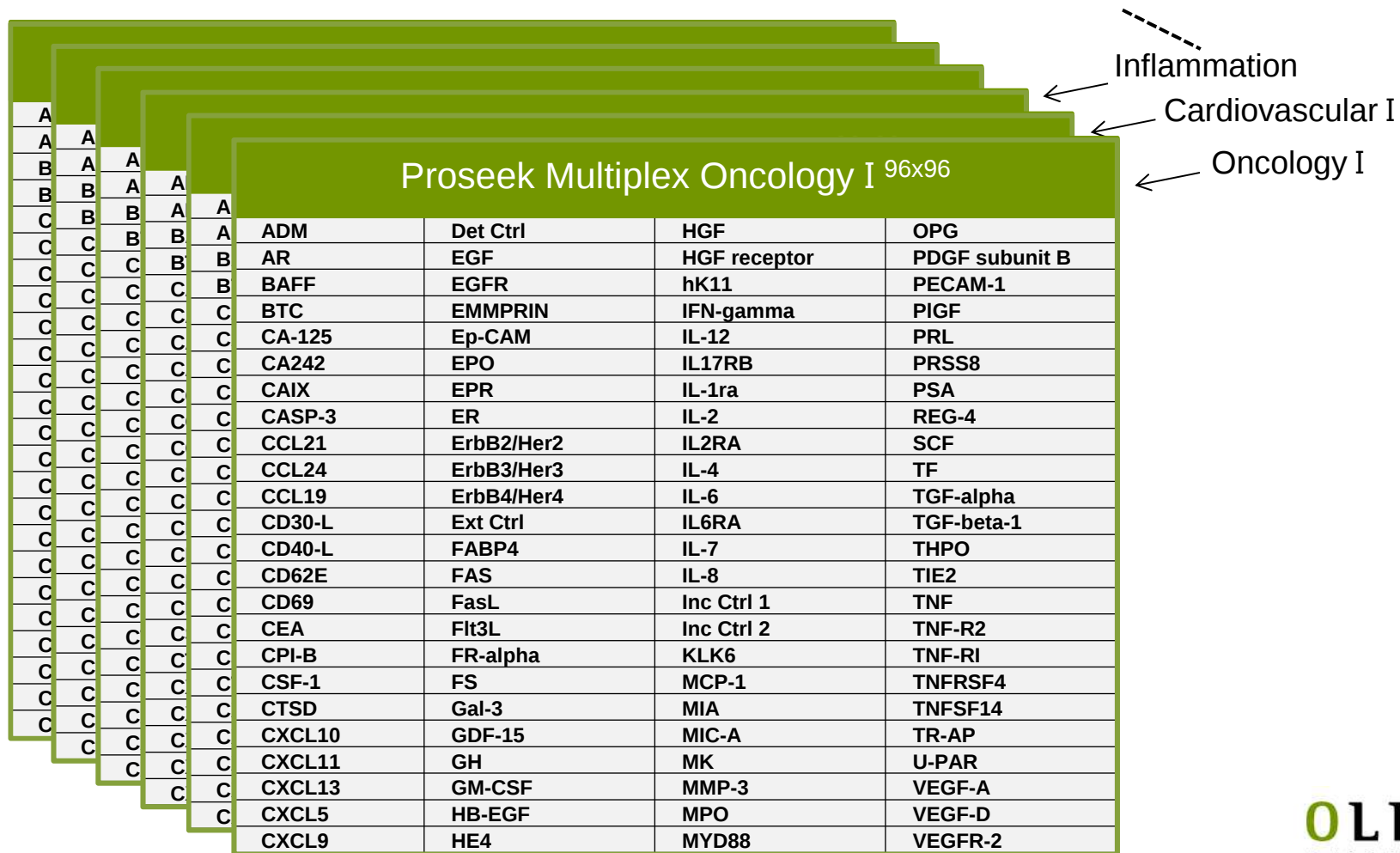

Proseek Multiplex Oncology I 96x96


ADM	Det Ctrl	HGF	OPG
AR	EGF	HGF receptor	PDGF subunit B
BAFF	EGFR	hK11	PECAM-1
BTC	EMMPRIN	IFN-gamma	PIGF
CA-125	Ep-CAM	IL-12	PRL
CA242	EPO	IL17RB	PRSS8
CAIX	EPR	IL-1ra	PSA
CASP-3	ER	IL-2	REG-4
CCL21	ErbB2/Her2	IL2RA	SCF
CCL24	ErbB3/Her3	IL-4	TF
CCL19	ErbB4/Her4	IL-6	TGF-alpha
CD30-L	Ext Ctrl	IL6RA	TGF-beta-1
CD40-L	FABP4	IL-7	THPO
CD62E	FAS	IL-8	TIE2
CD69	FasL	Inc Ctrl 1	TNF
CEA	Flt3L	Inc Ctrl 2	TNF-R2
CPI-B	FR-alpha	KLK6	TNF-RI
CSF-1	FS	MCP-1	TNFRSF4
CTSD	Gal-3	MIA	TNFSF14
CXCL10	GDF-15	MIC-A	TR-AP
CXCL11	GH	MK	U-PAR
CXCL13	GM-CSF	MMP-3	VEGF-A
CXCL5	HB-EGF	MPO	VEGF-D
CXCL9	HE4	MYD88	VEGFR-2

Inflammation
 Cardiovascular I
 Oncology I

Our first protein biomarker panel...of many yet to come


Proseek® Multiplex Oncology I ^{96x96}

Proseek Multiplex Oncology I ^{96x96}		
Adrenomedullin (ADM)	Fatty acid binding protein 4 adipocyte (FABP4)	Myeloperoxidase (MPO)
Amphiregulin (AR)	Fms-related tyrosine kinase 3 ligand (Flt3L)	Detection Control (Det Ctrl)
Angiopoietin-1 receptor (TIE2)	Folate receptor alpha (FR-alpha)	Osteopontin (OPN)
B-cell activating factor (BAFF)	Follistatin (FS)	Ovarian cancer-related tumor marker 125 (CA-125)
Betacellulin (BTC)	Galectin-3 (Gal-3)	Incubation Control (Inc Ctrl 2)
CA242 tumor marker (CA242)	Granulocyte-macrophage colony-stimulating factor (GM-CSF)	Placenta Growth Factor (PlGF)
Carbonic Anhydrase IX (CAIX)	Incubation Control (Inc Ctrl 1)	Platelet endothelial cell adhesion molecule (PECAM-1)
Carcinoembryonic antigen (CEA)	Growth Hormone (GH)	Platelet-derived growth factor subunit B (PDGF subunit B)
Caspase-3 (CASP-3)	Growth/differentiation factor 15 (GDF-15)	Prolactin (PRL)
Cathepsin D (CTSD)	Heparin-binding EGF-like growth factor (HB-EGF)	Prostate-specific antigen (PSA)
C-C motif chemokine 19 (CCL19)	Hepatocyte growth factor (HGF)	Prostate-specific antigen (PSA)
C-C motif chemokine 21 (CCL21)	Hepatocyte growth factor receptor (HGF receptor)	Receptor tyrosine-protein kinase ErbB-2 (ErbB2/Her2)
C-C motif chemokine 24 (CCL24)	Interferon gamma (IFN-gamma)	Receptor tyrosine-protein kinase ErbB-3 (ErbB3/Her3)
CD40 ligand (CD40-L)	Interleukin 1 receptor antagonist protein (IL-1ra)	Receptor tyrosine-protein kinase ErbB-4 (ErbB4/Her4)
C-X-C motif chemokine 10 (CXCL10)	Interleukin 12 (IL-12)	Regenerating islet-derived protein 4 (REG-4)
C-X-C motif chemokine 11 (CXCL11)	Interleukin 17 receptor B (IL17RB)	Stem cell factor (SCF)
C-X-C motif chemokine 13 (CXCL13)	Interleukin 2 (IL-2)	Tartrate-resistant acid phosphatase type 5 (TR-AP)
C-X-C motif chemokine 5 (CXCL5)	Interleukin 2 receptor subunit alpha (IL2RA)	Thrombopoietin (THPO)
C-X-C motif chemokine 9 (CXCL9)	Interleukin 4 (IL-4)	Tissue Factor (TF)
Cystatin B (CPI-B)	Interleukin 6 (IL-6)	Transforming growth factor alpha (TGF-alpha)
Early activation antigen CD69 (CD69)	Interleukin 6 receptor subunit alpha (IL6RA)	Transforming growth factor beta 1 (TGF-beta-1)
Epidermal growth factor (EGF)	Interleukin 7 (IL-7)	Tumor necrosis factor alpha (TNF)
Epidermal growth factor receptor (EGFR)	Interleukin 8 (IL-8)	Tumor necrosis factor ligand superfamily member 14 (TNFSF14)
Epidermal secretory protein E4 (HE4)	Kallikrein-11 (hK11)	Tumor necrosis factor ligand superfamily member 8 (CD30-L)
Eprex (EPR)	Kallikrein-6 (KLK6)	Tumor necrosis factor receptor 1 (TNF-R1)
Epithelial cell adhesion molecule (Ep-CAM)	Macrophage colony-stimulating factor 1 (CSF-1)	Tumor necrosis factor receptor 2 (TNF-R2)
Erythropoietin (EPO)	Matrix metalloproteinase-3 (MMP-3)	Tumor necrosis factor receptor superfamily member 4 (TNFSF4)
E-selectin (CD62E)	Melanoma-derived growth regulatory protein (MDR)	Tumor necrosis factor receptor superfamily member 6 (FAS)
Estrogen receptor (ER)	MHC class I polypeptide-related sequence A (MIC-A)	Urokinase plasminogen activator surface receptor (U-PA)
Extension Control (Ext Ctrl)	Midkine (MK)	Vascular endothelial growth factor A (VEGF-A)
Extracellular matrix metalloproteinase inducer (EMMPRN)	Monocyte chemoattractant protein-1 (MCP-1)	Vascular endothelial growth factor D (VEGF-D)
Fas antigen ligand (FasL)	Myeloid differentiation primary response protein MyD88 (MYD88)	Vascular endothelial growth factor receptor 2 (VEGFR-2)

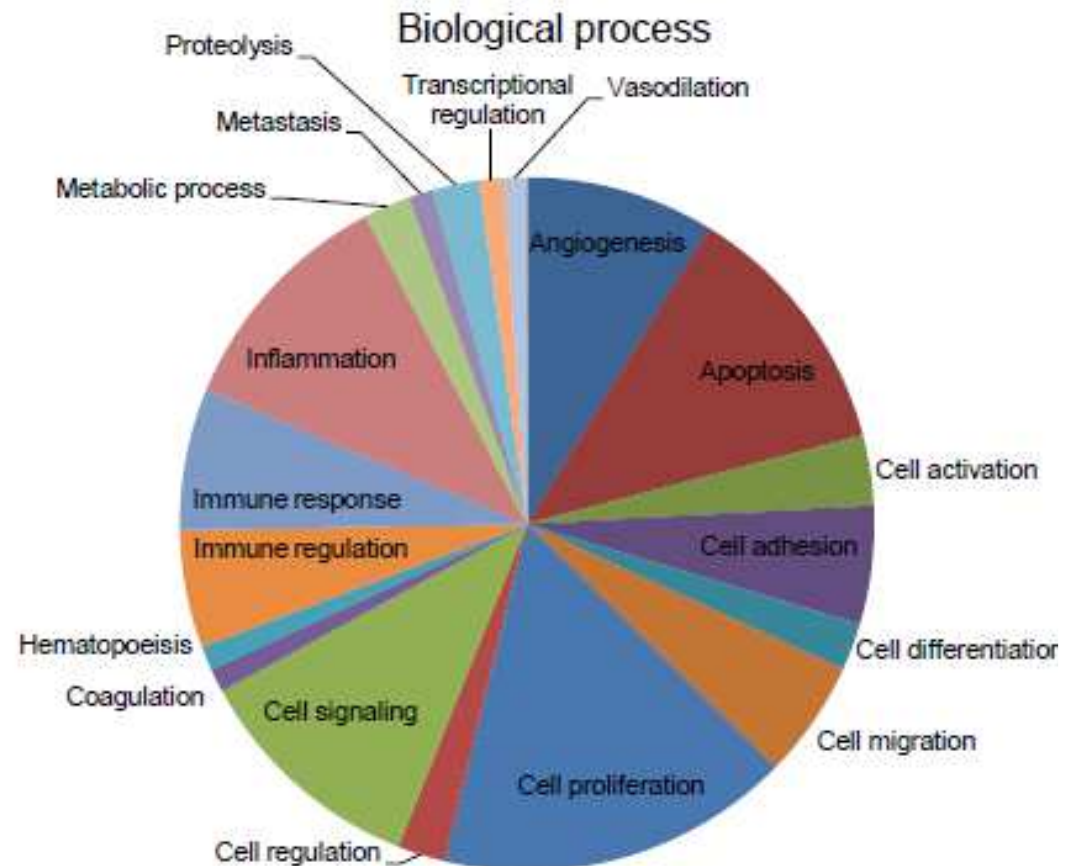


Real life example 1

- ▶ Markers for early detection of CRC
- ▶ Study performed with Proseek Multiplex Oncology I 96x96
- ▶ Thorsen et al submitted to Molecular and Cellular Proteomics

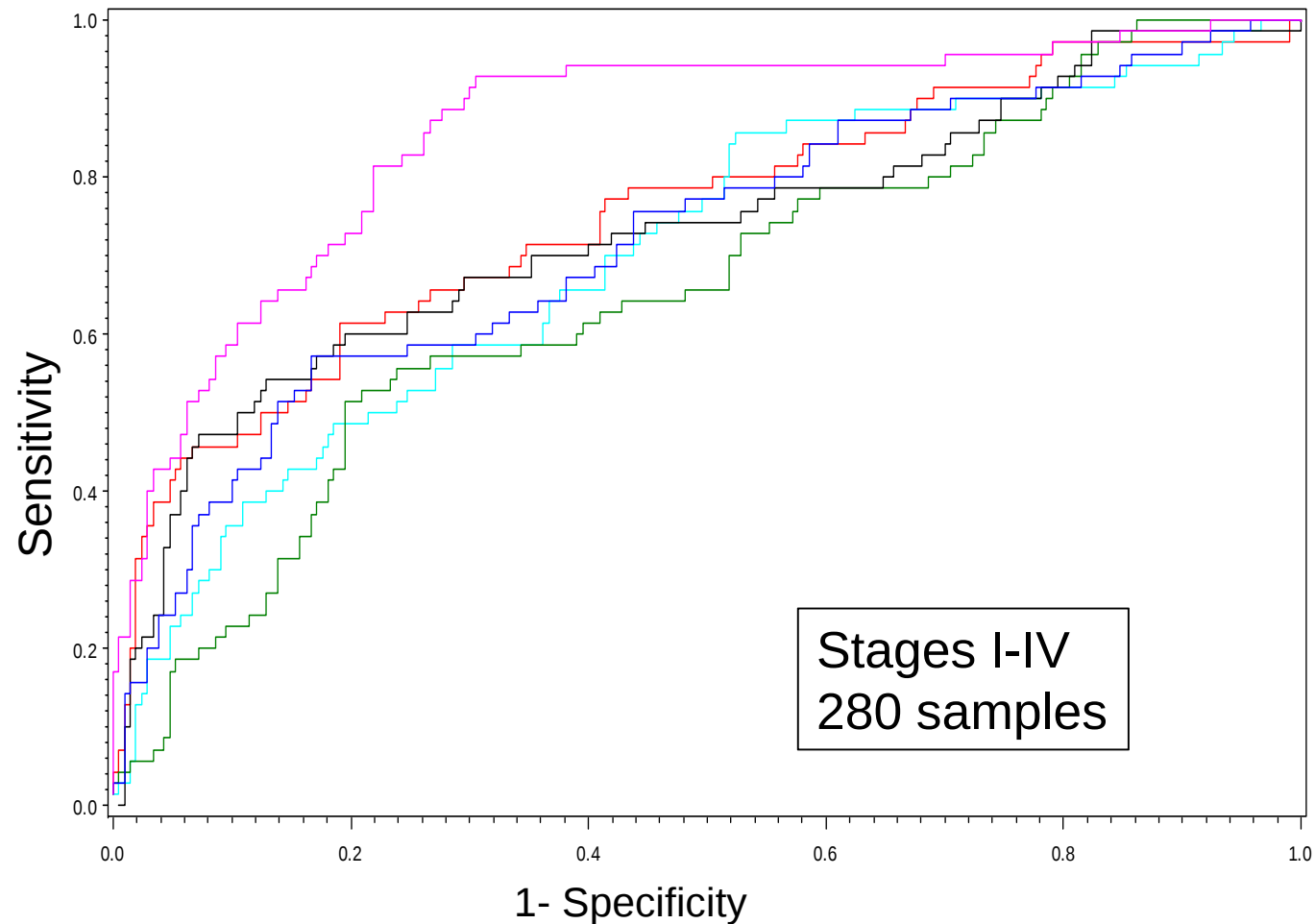





Real life example 2

Comparing groups of healthy and CVD patients

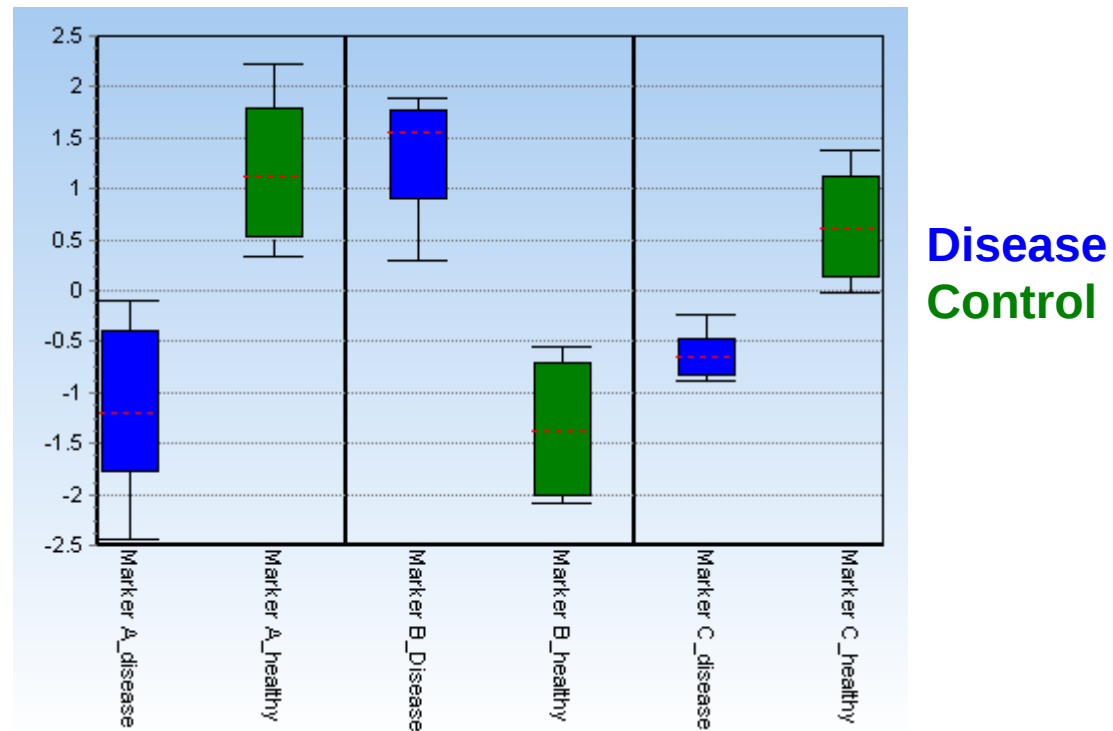



Real Life example 3

Comparing protein profiles over time in brain microdialysis fluid

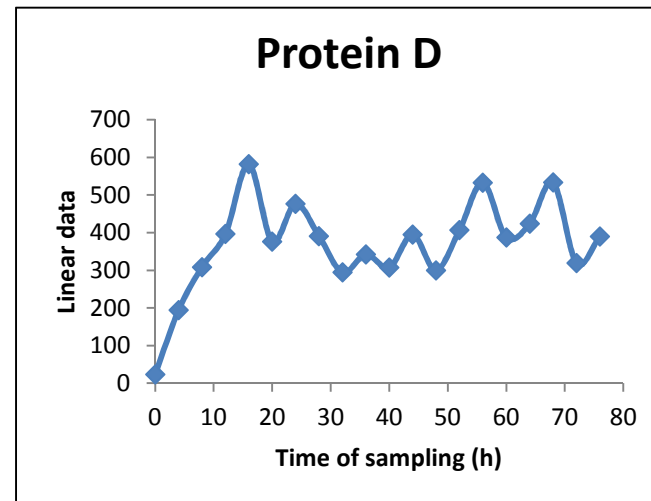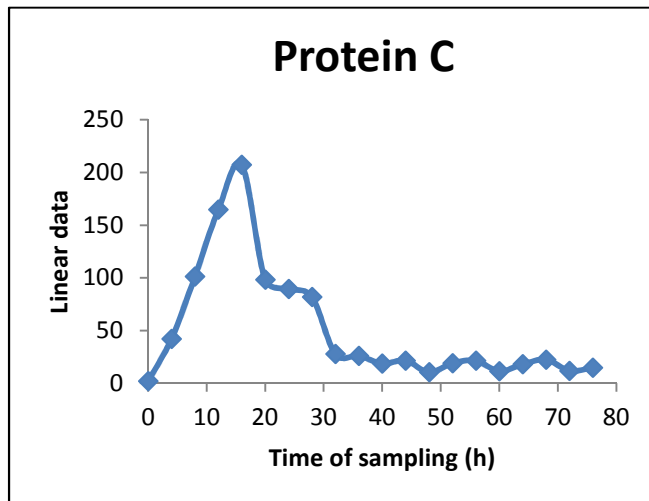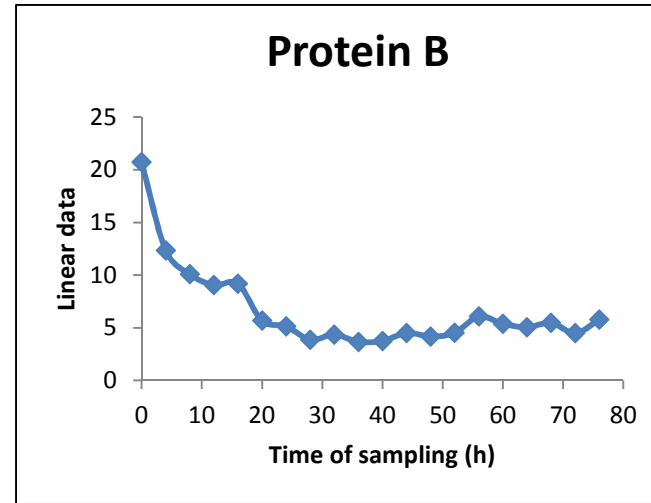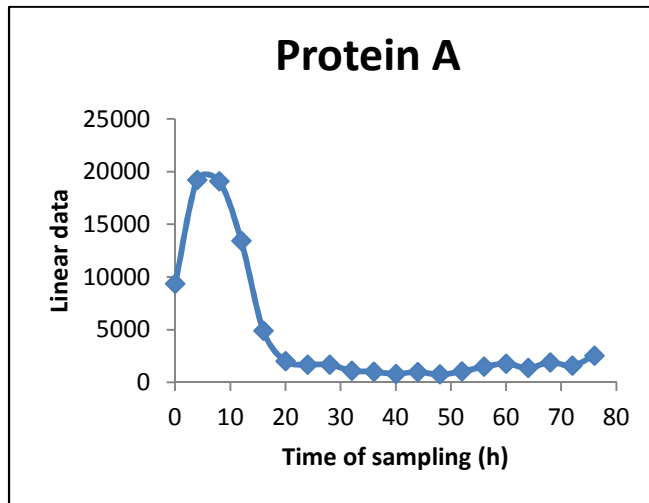

Real life example 4

Drug response biomarkers in Chronic Myelogenous Leukemia visualized with Principle Component Analysis and box plot

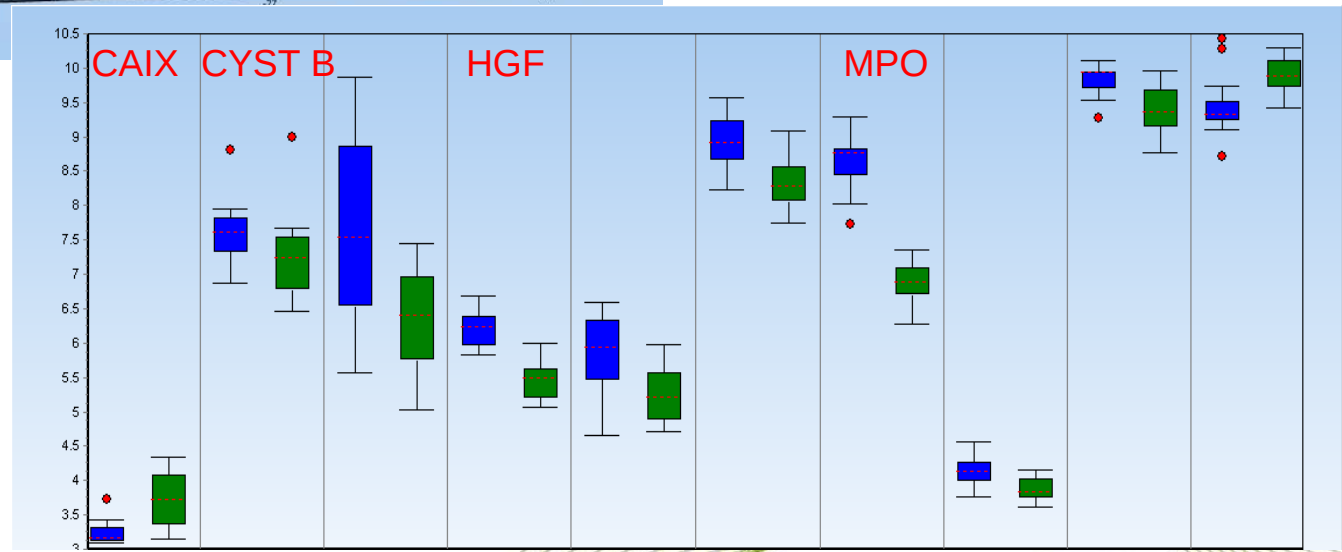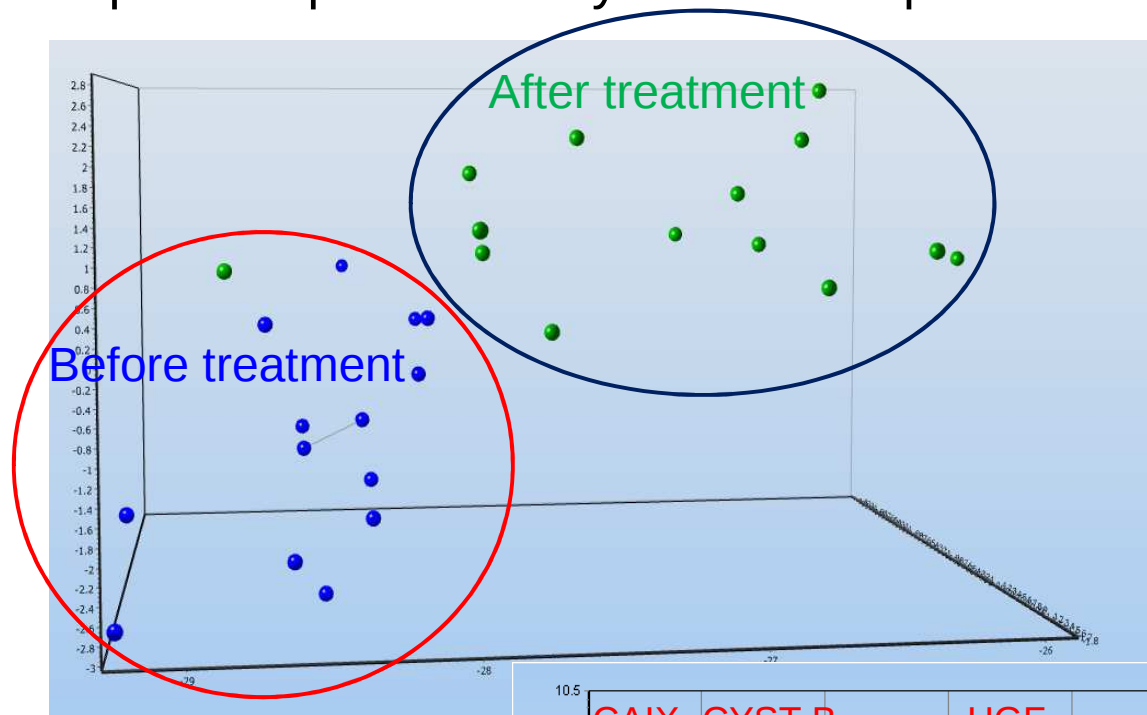

Data analysis software, customized for Proseek Multiplex

- ▶ Direct import of data from Biomark
- ▶ Has QC parameters and several statistical analysis methods
- ▶ <http://multid.se/OlinkWizardUserGuide.pdf>
- *The only product specific analysis tool for multivariate statistics in immunoassays*

Olink Wizard for GenEx

USER GUIDE

Version 1.0 (June 2013)



Paired t-tests identifies putative biomarkers

(A) vs (B)	Normality Tes...	Fold change	Difference (A-...	P-Value
MPO	Passed	3.4058	1.76799	1E-8
HGF	Passed	1.64382	0.71705	6.3608761975...
IL6RA	Passed	1.49349	0.57869	9.5423780435...
E-selectin	Passed	2.70588	1.4361	3.5746934282...
VEGF-D	Passed	-1.38577	-0.47069	0.0001312826...
U-PAR	Passed	1.35344	0.43663	0.0001983101...
IL-1ra	Passed	1.49963	0.58461	0.0002112718...
Cystatin B	Passed	1.27812	0.35403	0.0004590326...
TNFSF14	Passed	1.19635	0.25864	0.0005459262...
CAIX	Passed	-1.41649	-0.50232	0.0005661618...
TNF-RI	Passed	1.25934	0.33266	0.0006719713...
PIGF	Passed	1.288	0.36513	0.0012267408...
TNF-R2	Passed	1.23019	0.29888	0.0022588399...
IL2RA	Passed	1.15688	0.21023	0.0027132323...
Stem cell factor	Passed	-1.23166	-0.3006	0.0027877186...
FR-alpha	Passed	-1.08439	-0.11689	0.0033619099...
IL-12	Passed	-1.74466	-0.80295	0.0041622586...
TGF-alpha	Passed	1.32551	0.40654	0.0047751327...
ErbB4/Her4	Passed	-1.18967	-0.25057	0.0048166776...
PECAM-1	Passed	1.1673	0.22318	0.0073897061...
Kallikrein-11	Passed	-1.19558	-0.25771	0.0115430763...
IL-8	Passed	-1.58315	-0.6628	0.0121299853...
IL-10	Passed	1.12584	0.171	0.0150805018...

Bonferroni multiple correctio


$p < 0.00056$

$p > 0.00056$



Proseek Multiplex Technology


The technology is based on Proximity Extension Assay (PEA)

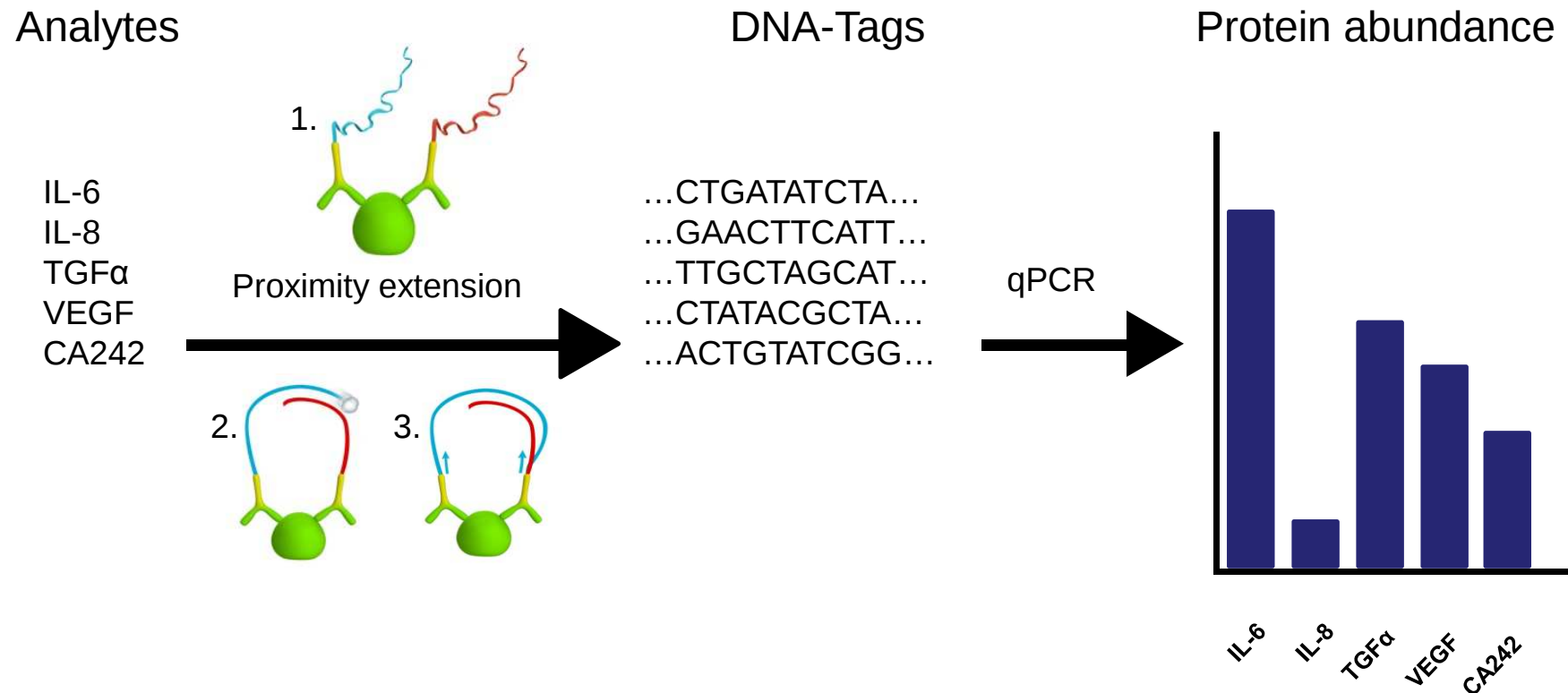


"A cDNA synthesis for proteins"

Multiplexing often means specificity problems

Problem: Multiplex sandwich immunoassays → cross reactivity

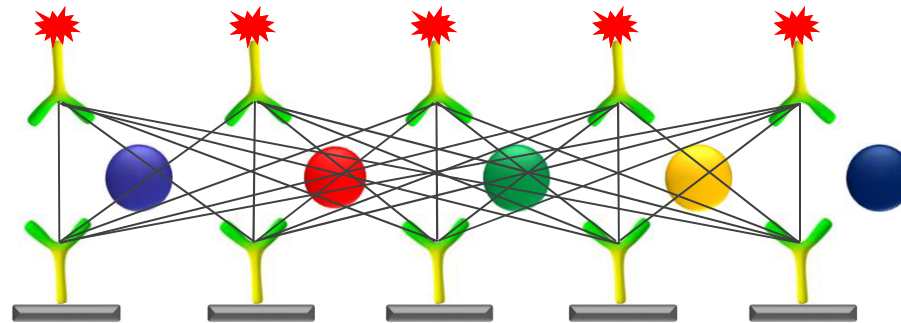

Solution: Only matched pairs detected through DNA-barcoding

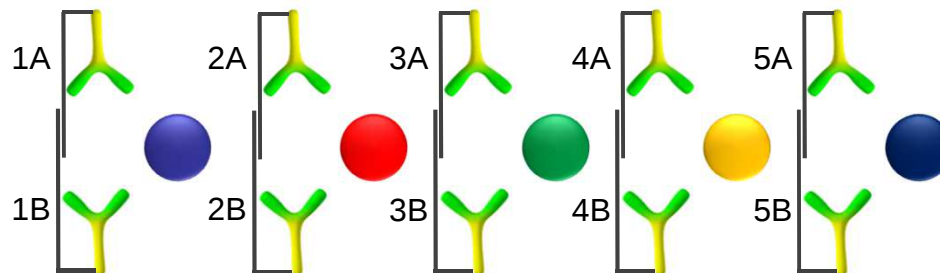

It works for different biological materials

- and only in 1 µl

- ▶ Biobanked material
 - ▶ Precious samples
 - ▶ Limited sample volume
- Serum and plasma
 - Tissue lysate
 - Cell lysate
 - Fine Needle Biopsy
 - Microdialysis fluid
 - Cell culture media
 - Serum from xenograft mice


"I really think that one of the best advantages of your platform is the low amount of sample required. This is a huge positive since many of our samples are quite precious."

Sharon Pitteri, Stanford, Canary foundation

How does this sound?


- ▶ True scalability
- ▶ Small sample consumption
- ▶ No cross-reactive events detected
- ▶ Many different sample types

.....there is more

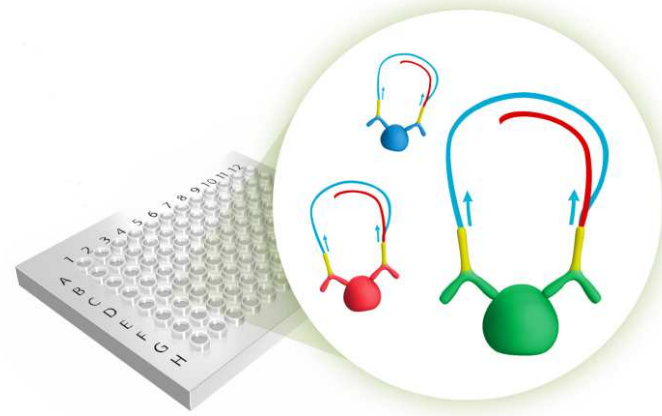


It is not even complicated


PREPARATION

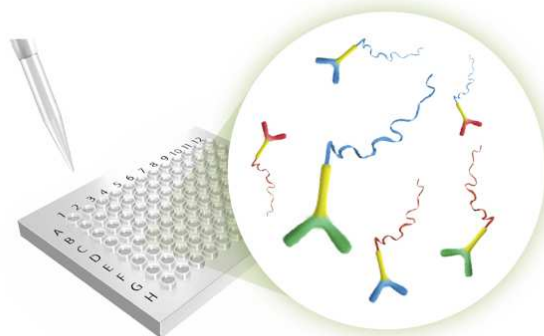

INCUBATION


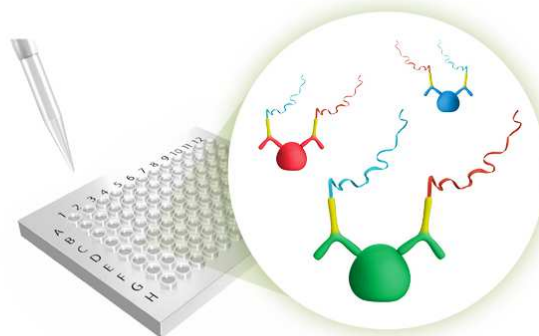


EXTENSION

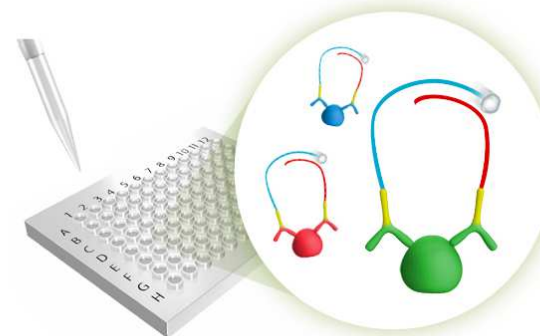

DETECTION


SAMPLE ANALYSIS

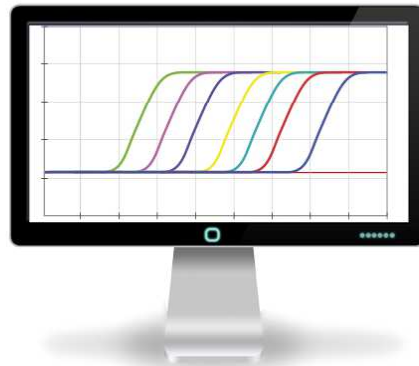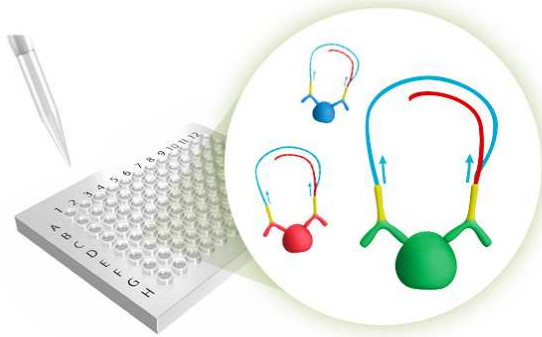

Hands-on-time: less than 3 hours



Four internal controls are included in every assay

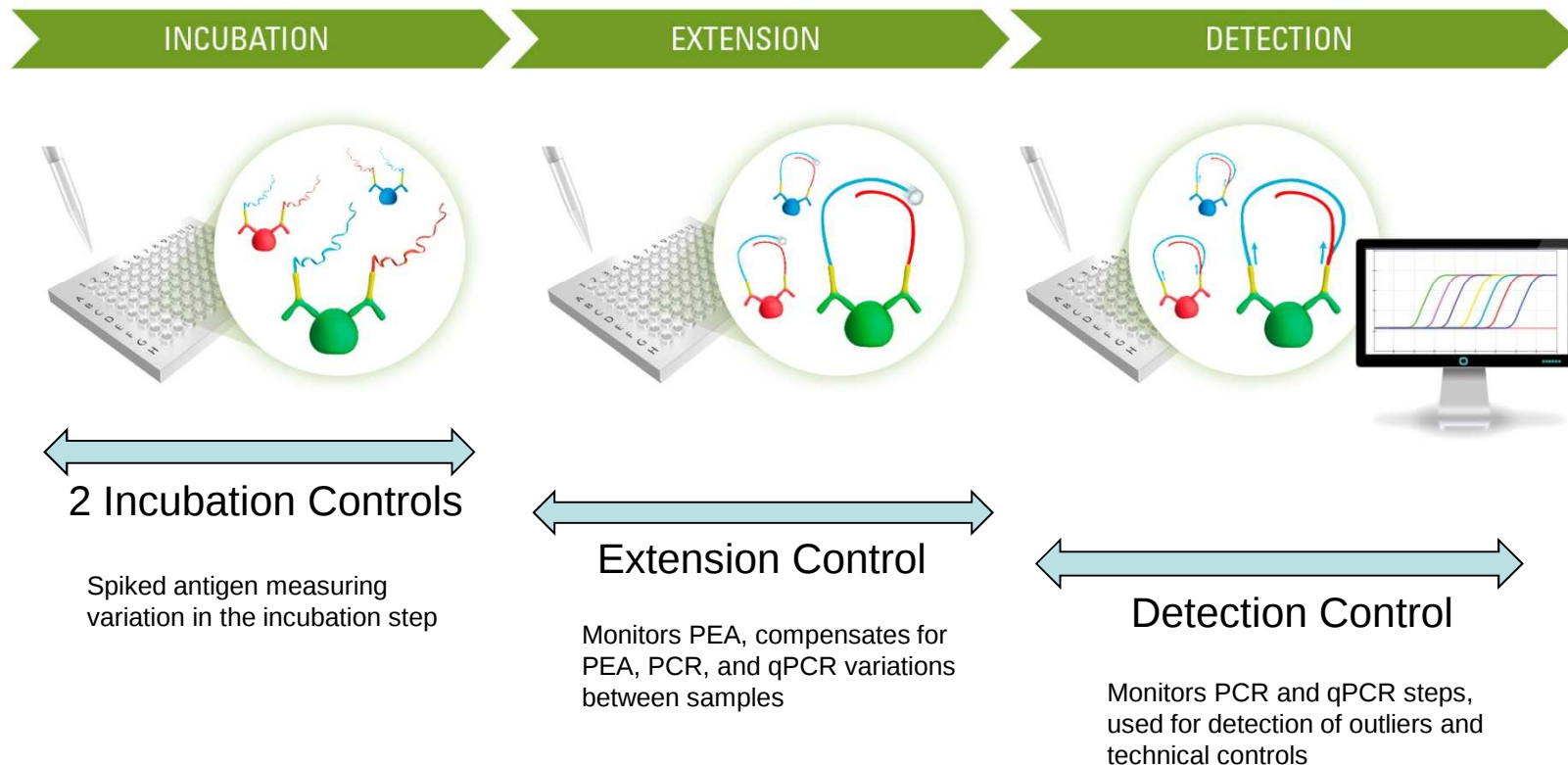


We use high throughput real time PCR





- ▶ Microfluidic system – Fluidigm Biomark system
- ▶ 96.96 Dynamic Array™ chip measuring 96x96 reactions
- ▶ 9,216 data points per run in 7 nL reaction chambers

Typical data readout

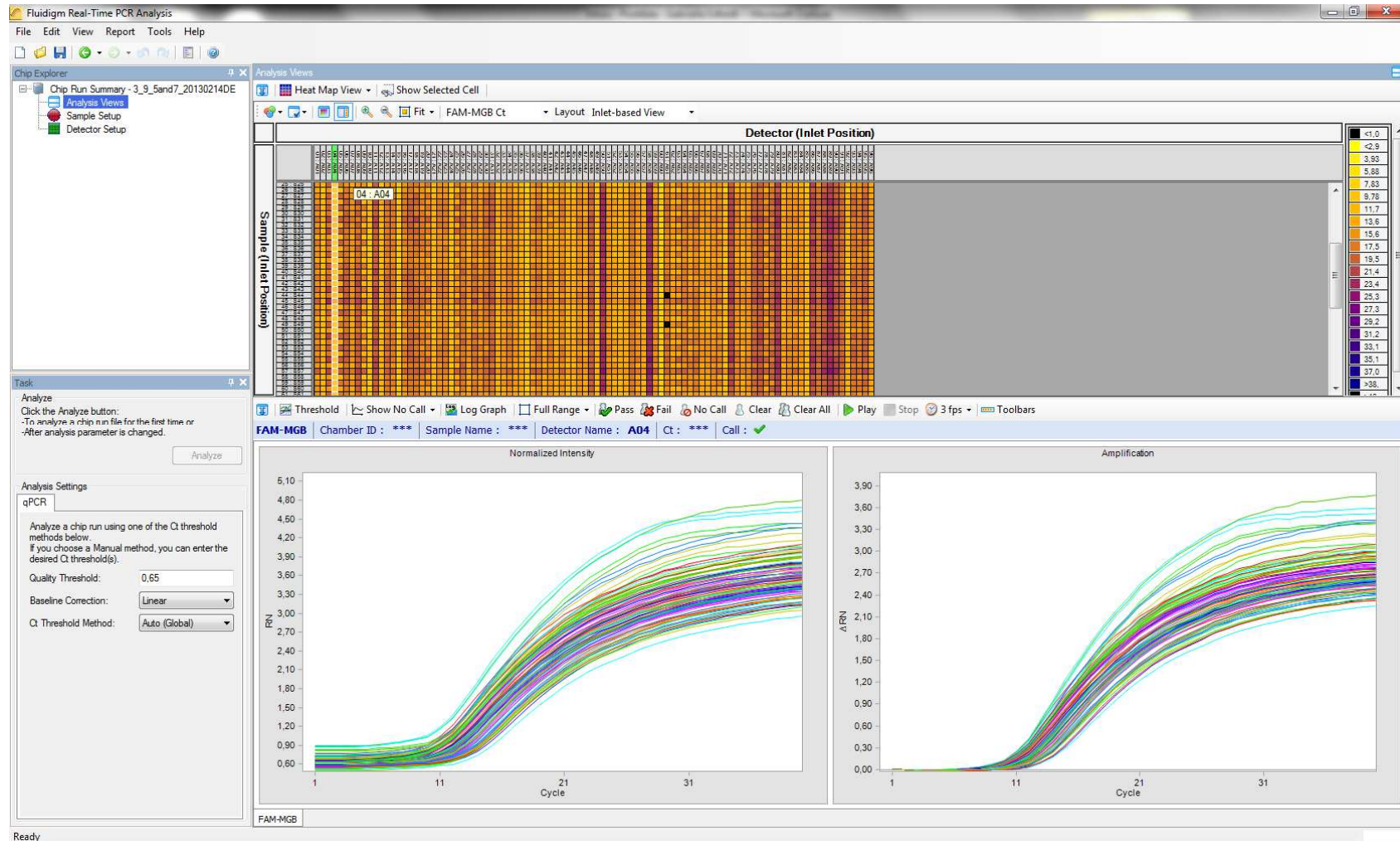

Olink Wizard in GenEx helps to analyze data

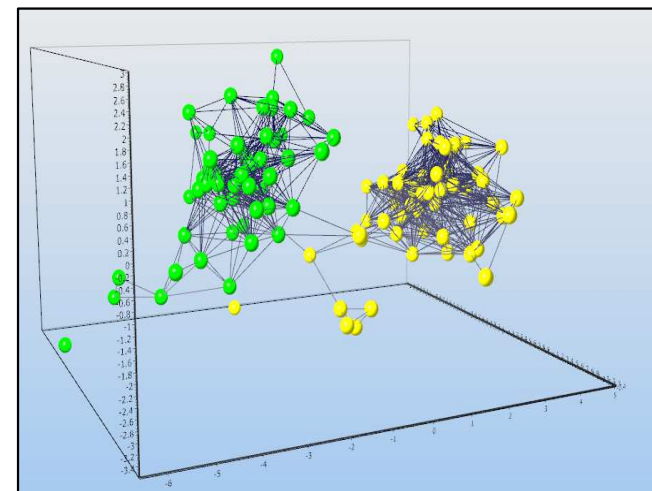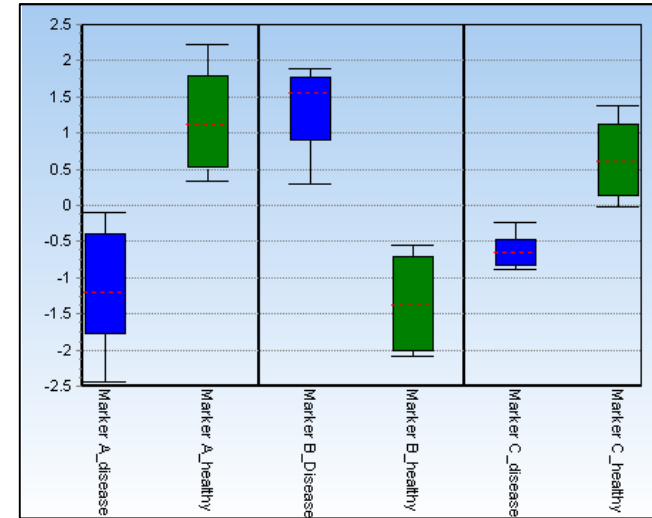

K

BIOSCIENCE


The product

- ▶ Proseek Multiplex Probe Kit 96x96
- ▶ Proseek Multiplex Detection Kit 96x96
- ▶ Proseek Multiplex Controls



www.olink.com

Coming next.....

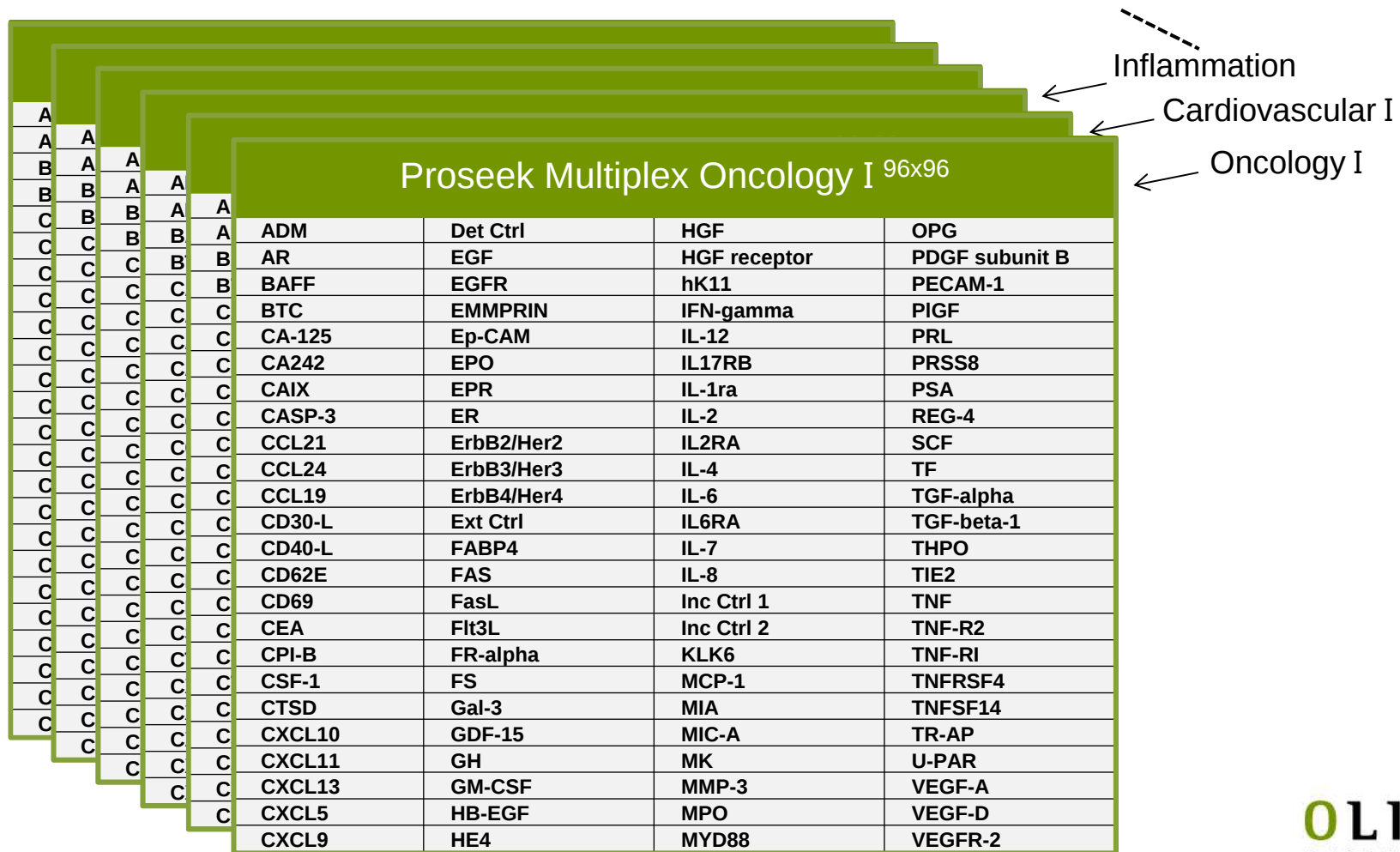

Proseek Multiplex Oncology I 96x96


ADM	Det Ctrl	HGF	OPG
AR	EGF	HGF receptor	PDGF subunit B
BAFF	EGFR	hK11	PECAM-1
BTC	EMMPRIN	IFN-gamma	PIGF
CA-125	Ep-CAM	IL-12	PRL
CA242	EPO	IL17RB	PRSS8
CAIX	EPR	IL-1ra	PSA
CASP-3	ER	IL-2	REG-4
CCL21	ErbB2/Her2	IL2RA	SCF
CCL24	ErbB3/Her3	IL-4	TF
CCL19	ErbB4/Her4	IL-6	TGF-alpha
CD30-L	Ext Ctrl	IL6RA	TGF-beta-1
CD40-L	FABP4	IL-7	THPO
CD62E	FAS	IL-8	TIE2
CD69	FasL	Inc Ctrl 1	TNF
CEA	Flt3L	Inc Ctrl 2	TNF-R2
CPI-B	FR-alpha	KLK6	TNF-RI
CSF-1	FS	MCP-1	TNFRSF4
CTSD	Gal-3	MIA	TNFSF14
CXCL10	GDF-15	MIC-A	TR-AP
CXCL11	GH	MK	U-PAR
CXCL13	GM-CSF	MMP-3	VEGF-A
CXCL5	HB-EGF	MPO	VEGF-D
CXCL9	HE4	MYD88	VEGFR-2

Inflammation
Cardiovascular I
Oncology I

Cardio 96x96 (coming in October)

Adrenomodulin	DKK-1	IL-10	MMP-3	Thrombomodulin
AgRP	ECP	IL-17	MMP-7	Tie-2
Aminopeptidase P2	EGF	IL-1alpha	MMP-9	TIM-1
bFGF	Endoglin	IL-1beta	MPO	Tissue Factor
bNGF	EPO	IL-1RA	Myoglobin	TNF-R1
BNP	E-Selectin/CD62E	IL-2	NT-pro-BNP	TNF-R2
CA125	ESM-1	IL-25	Osteoprotegerin	tPA
Cathepsin D	FABP3	IL-4	PaPPA	TRAIL
Cathepsin L	FABP4	IL-6	PDGF-BB	TRANCE
Cathepsin S	Fas	IL-6R	Pecam-1/CD31	U-PAR
CCL3	FGF-23	IL-8	PLA2G7	VEGF
CCL4	Follistatin	Kallikrein 11	PIGF	VEGF D
CD40	F-Spondin	Kallikrein-6	Prolactin	YKL-40
CD40L	Galanin	LIGHT	PSGL-1	
CX3CL1	Galectin-3	MCP-1	PTX3	
CXCL-16	GDF-15	M-CSF	RAGE	
Cystatin B	GH	MMP-1	Renin	
DKK-1	HB-EGF	MMP-10	Resistin	
	HGF	MMP-12	SCF	
	HSP27		ST2	

IL-18

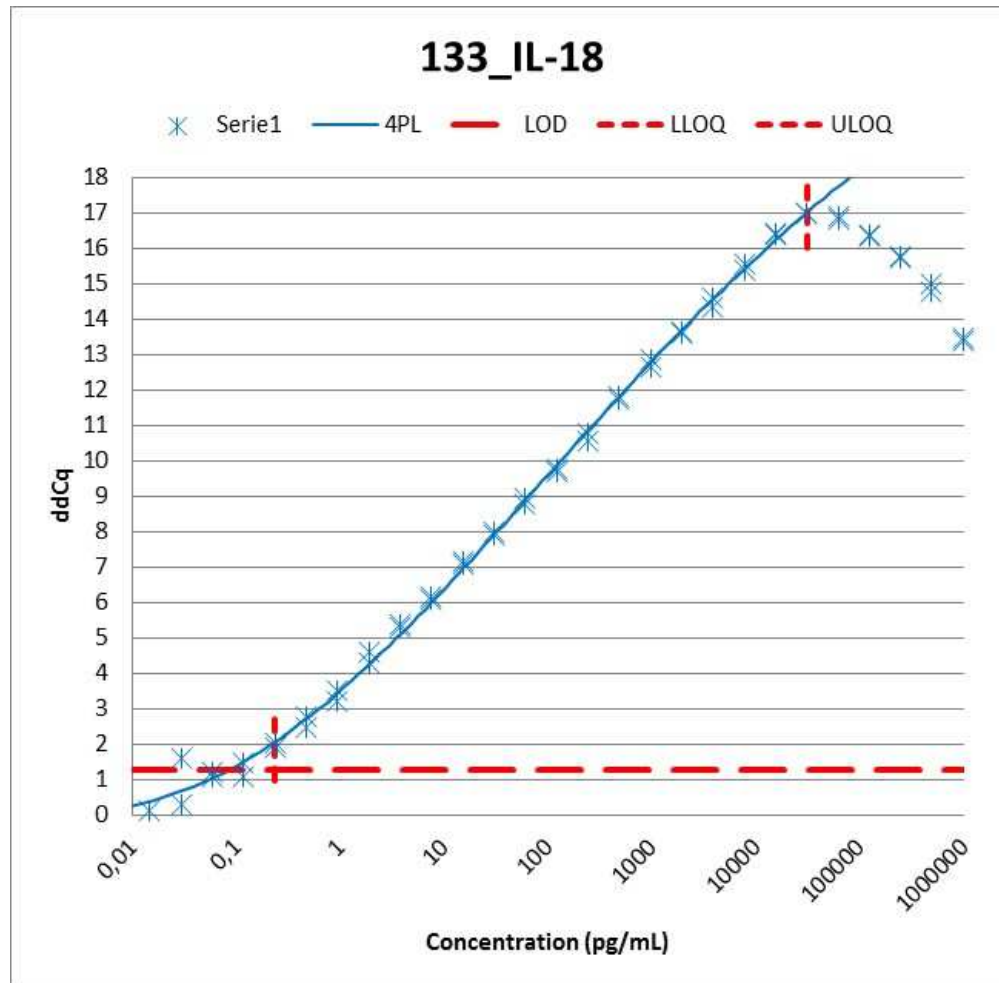


Recovery:91%

**No interference from:
Lipid, Haemolysate, Bilirubin**

LOD	0,12	pg/mL
LLOQ	0,24	pg/mL
ULOQ	31250	pg/mL
Range	5,12	log10

IL-4

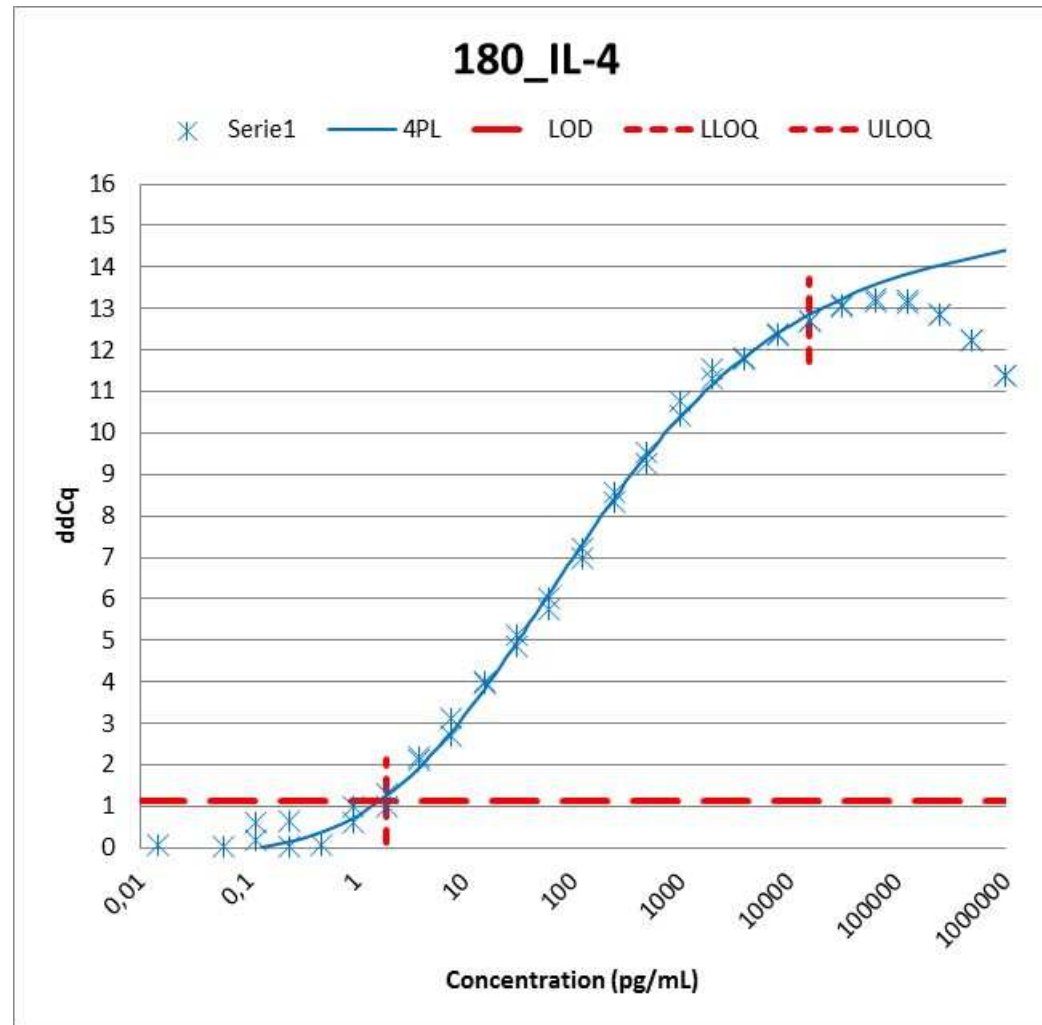


Recovery:105%

**No interference from:
Lipid, Haemolysate, Bilirubin**

LOD	1,91	pg/mL
LLOQ	1,91	pg/mL
ULOQ	15630	pg/mL
Range	3,91	log10

IL-6

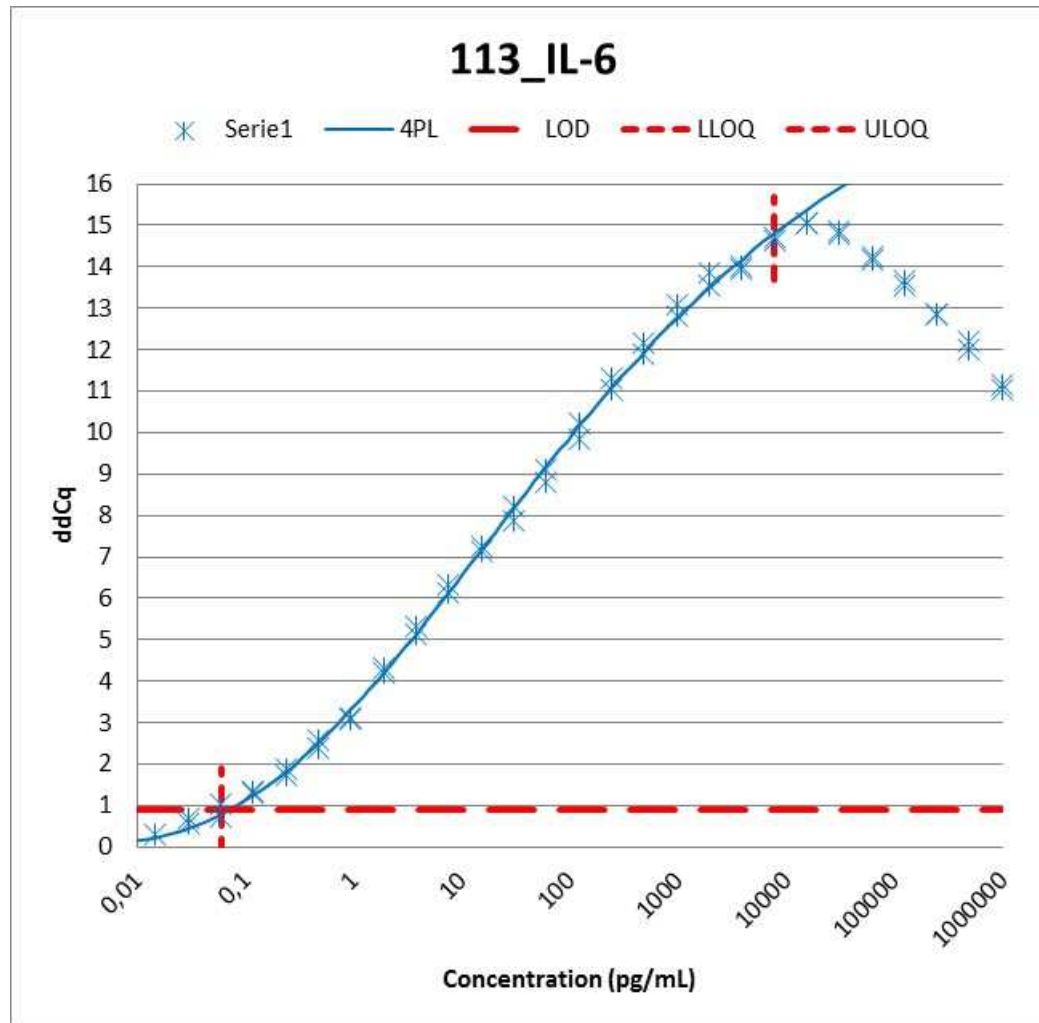


Recovery:108%

**No interference from:
Lipid, Haemolysate, Bilirubin**

LOD	0,06	pg/mL
LLOQ	0,06	pg/mL
ULOQ	7810	pg/mL
Range	5,12	log10

IL-6

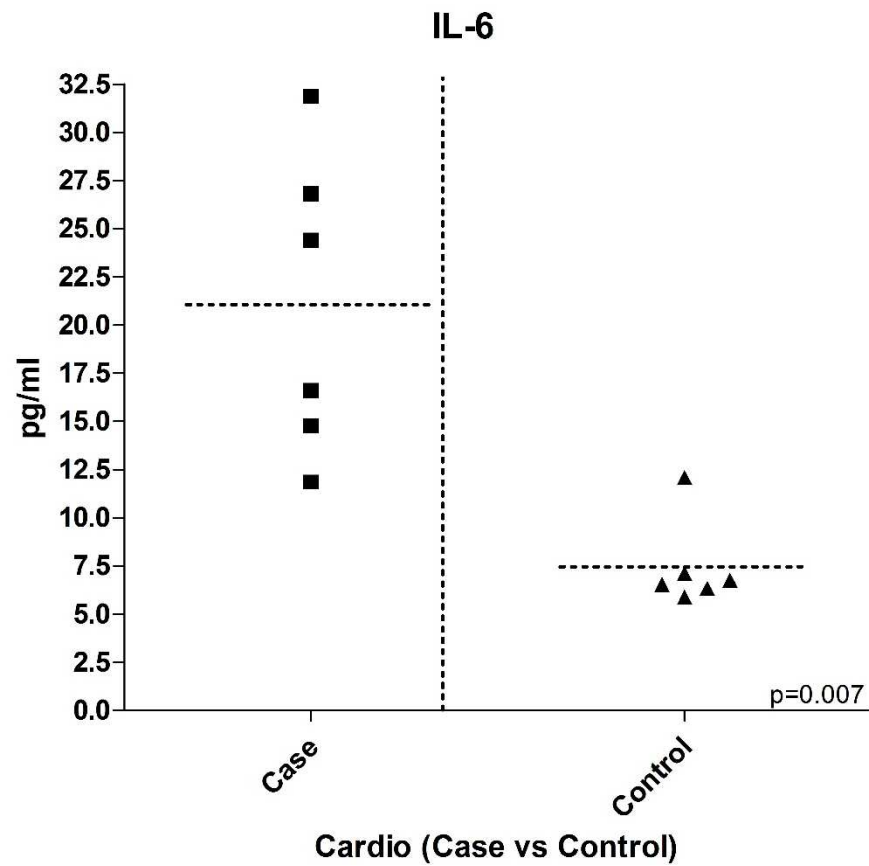


Recovery:108%

**No interference from:
Lipid, Haemolysate, Bilirubin**

LOD	0,06	pg/mL
LLOQ	0,06	pg/mL
ULOQ	7810	pg/mL
Range	5,12	log10

EGF

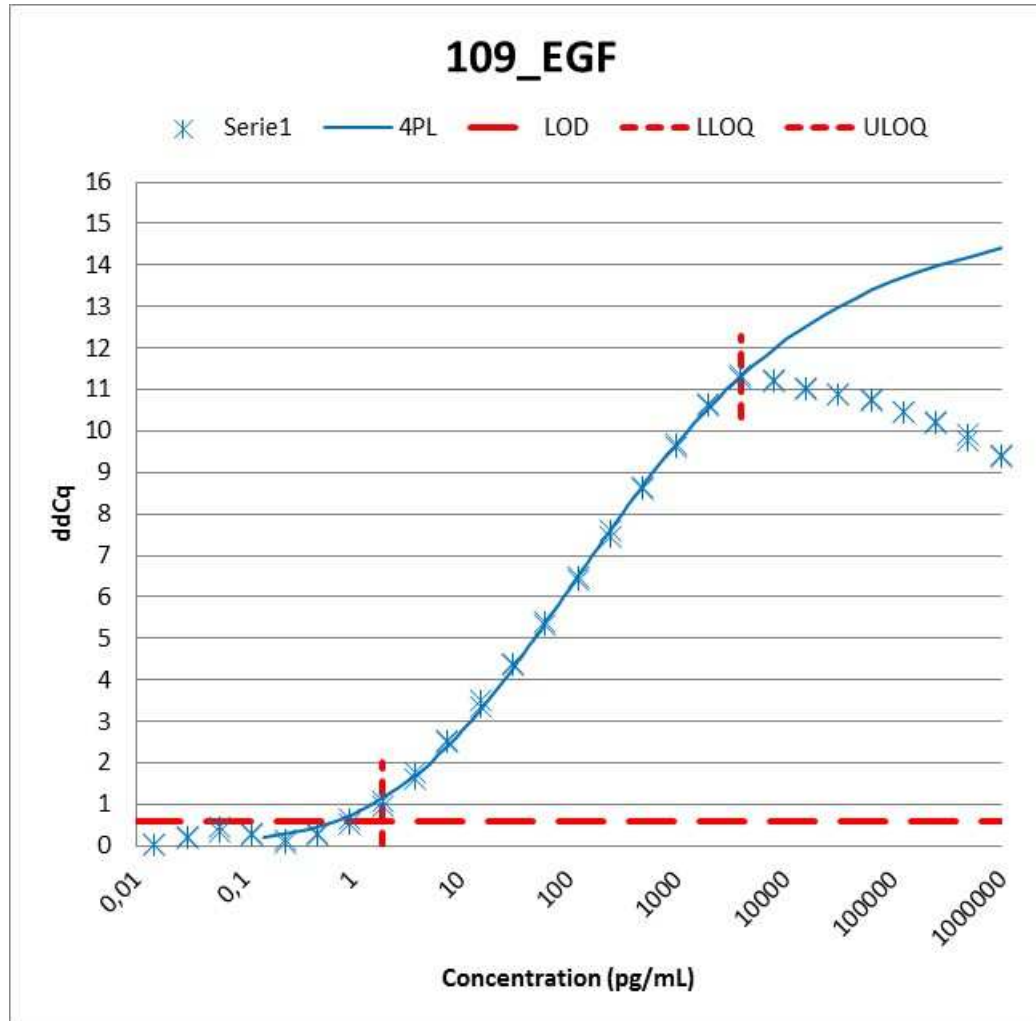


Recovery:100%

No interference from:
Lipid, Haemolysate, Bilirubin

LOD	0,95 pg/mL
LLOQ	1,91 pg/mL
ULOQ	3910 pg/mL
Range	3,31 log10

EGF

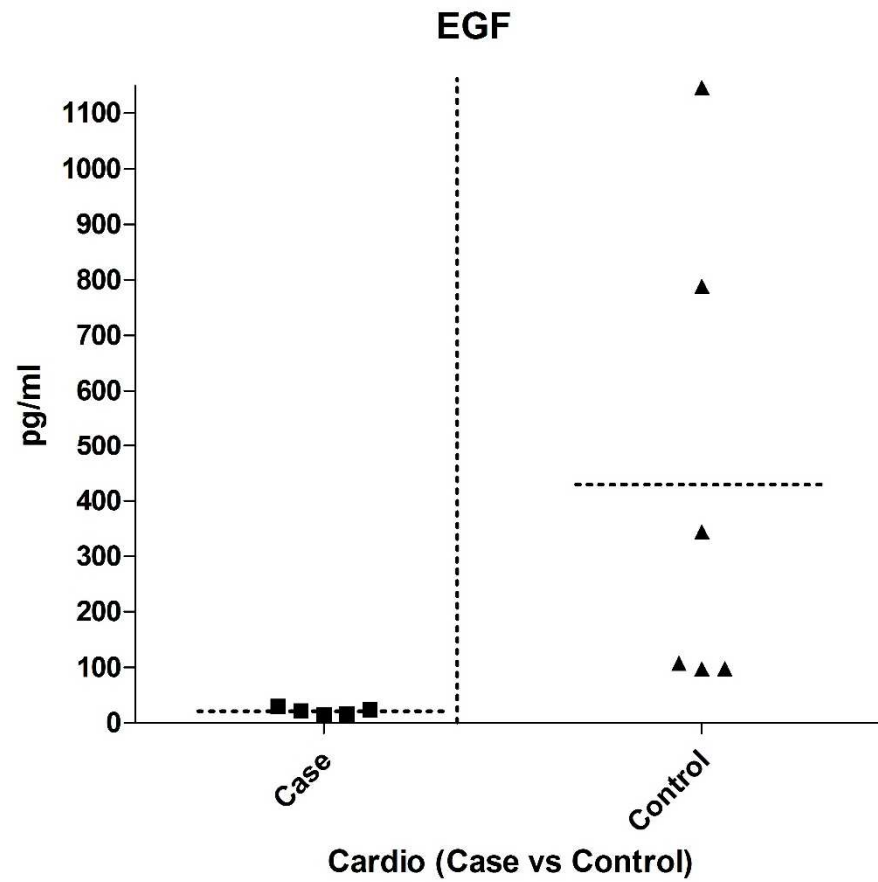


Recovery:100%

**No interference from:
Lipid, Haemolysate, Bilirubin**

LOD	0,95 pg/mL
LLOQ	1,91 pg/mL
ULOQ	3910 pg/mL
Range	3,31 log10

CXCL6

168_CXCL6

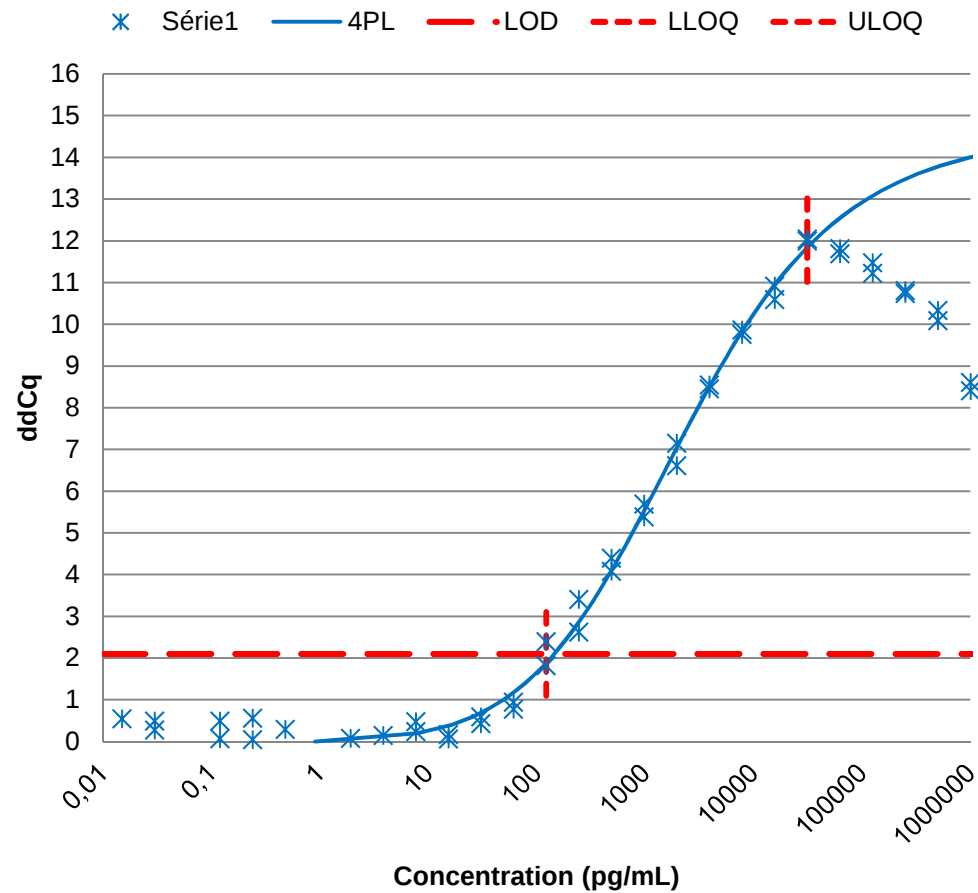


Recovery:95%

No interference from:
Lipid, Haemolysate, Bilirubin

LOD	122,07	pg/mL
LLOQ	122,07	pg/mL
ULOQ	31250	pg/mL
Range	2,41	log10

CXCL6

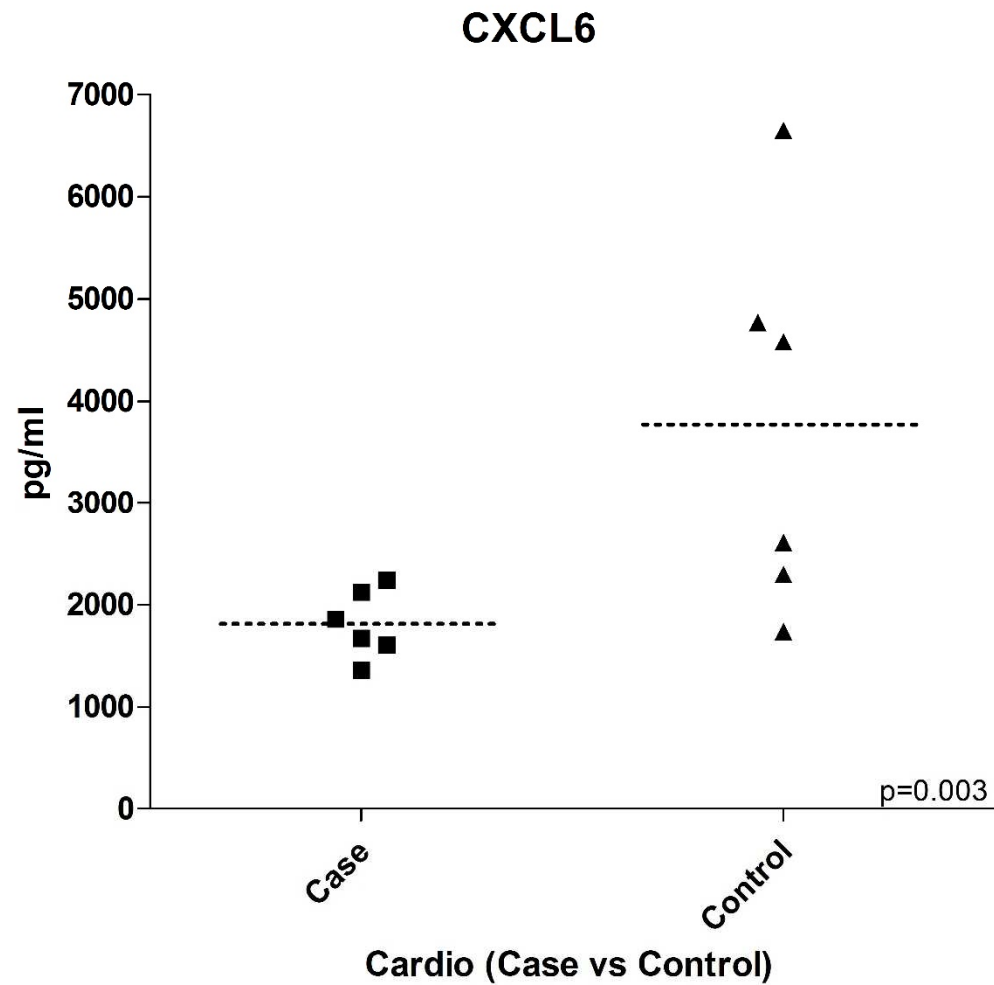


Recovery:95%

**No interference from:
Lipid, Haemolysate, Bilirubin**

LOD	122,07	pg/mL
LLOQ	122,07	pg/mL
ULOQ	31250	pg/mL
Range	2,41	log10

MMP-10

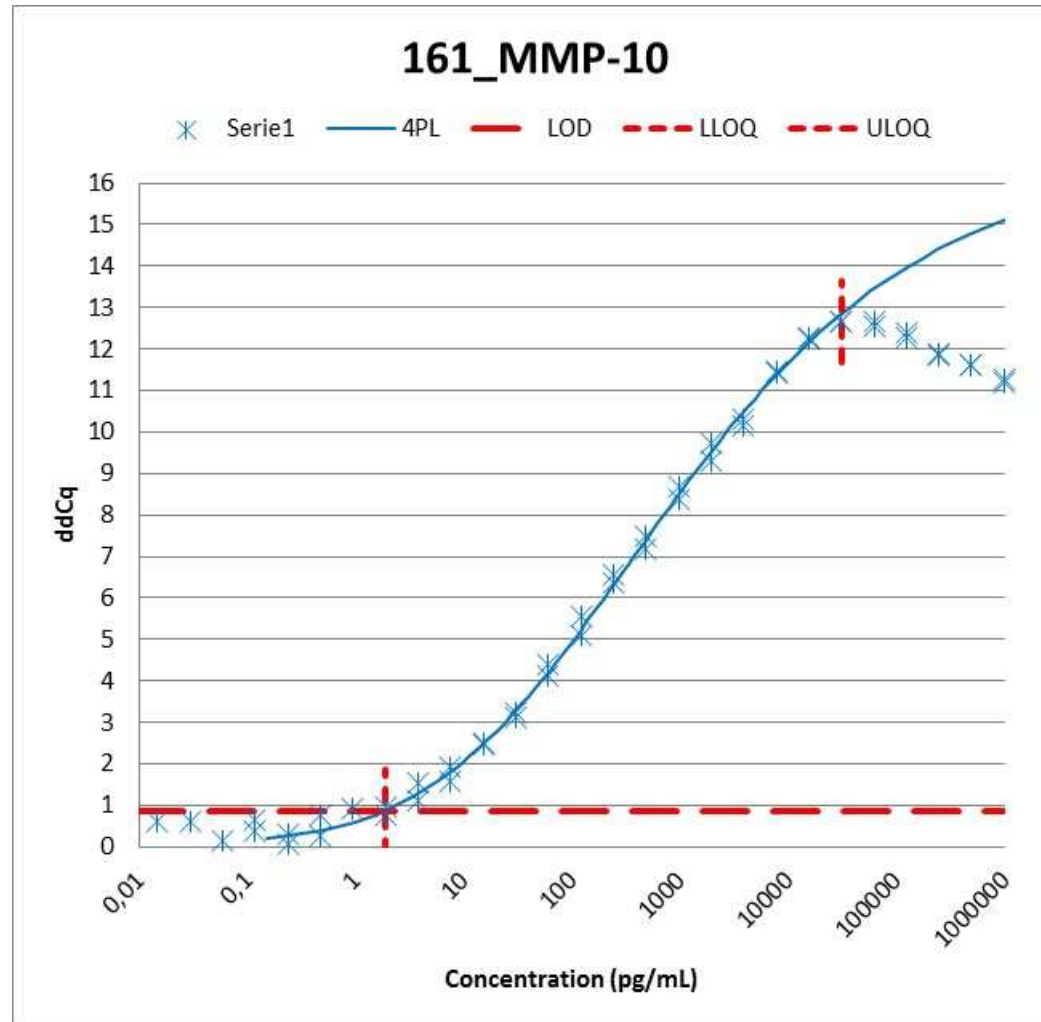

Recovery:100%

**No interference from:
Lipid, Haemolysate, Bilirubin**

LOD	1,91	pg/mL
LLOQ	1,91	pg/mL
ULOQ	31250	pg/mL
Range	4,21	log10

Who we are



- ▶ Founded in 2004
- ▶ Headquartered in Uppsala, Sweden
- ▶ Privately held
- ▶ Serving the top 10 global pharma
- ▶ Propriety immunoassay platform
- ▶ Proseek[®] & Duolink[®] Product Lines
(> 1000 users worldwide, > 35 patents and > 500 scientific publications)

www.olink.com

Proseek Multiplex -Scalable Immunoassays

Available at:

Olink Bioscience

Contact for Europe

Mats Bergström

mats.bergstrom@olink.com

Phone : +46 18 444 39 70

E-mail : order@olink.com



Performance

<http://www.olinK.com/products/proseek-multiplex/downloads/data-packages>

Performance – Sample material

- ▶ **Complex biological samples**

- Serum
- Plasma (EDTA, Citrate, Heparin)


- ▶ **Additional biological samples tested as part of proof of principle studies**

- Fine needle aspiration biopsies
- Micro dialysis fluid
- Tissue lysate
- Cell culture media
- Cell lysate
- Serum from xenograft mice


Biological material is complex

For example blood:

- ▶ Over 100 000 different proteins
- ▶ 22 most abundant proteins = 99% of the content

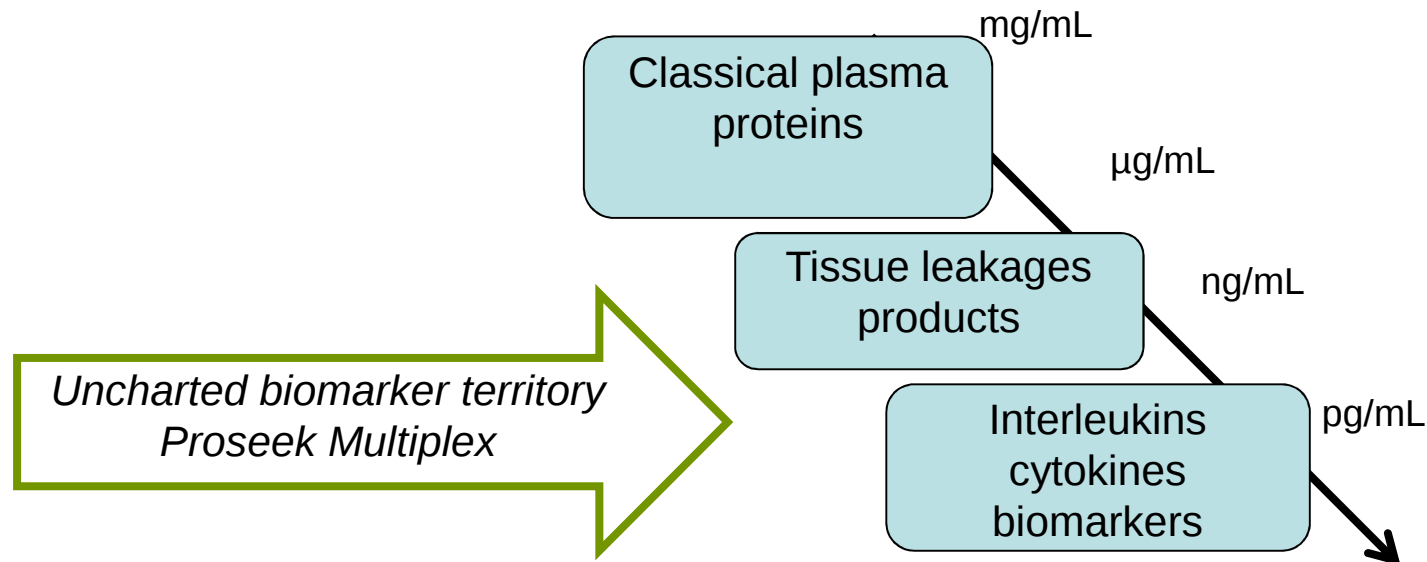

Performance – Dynamic range

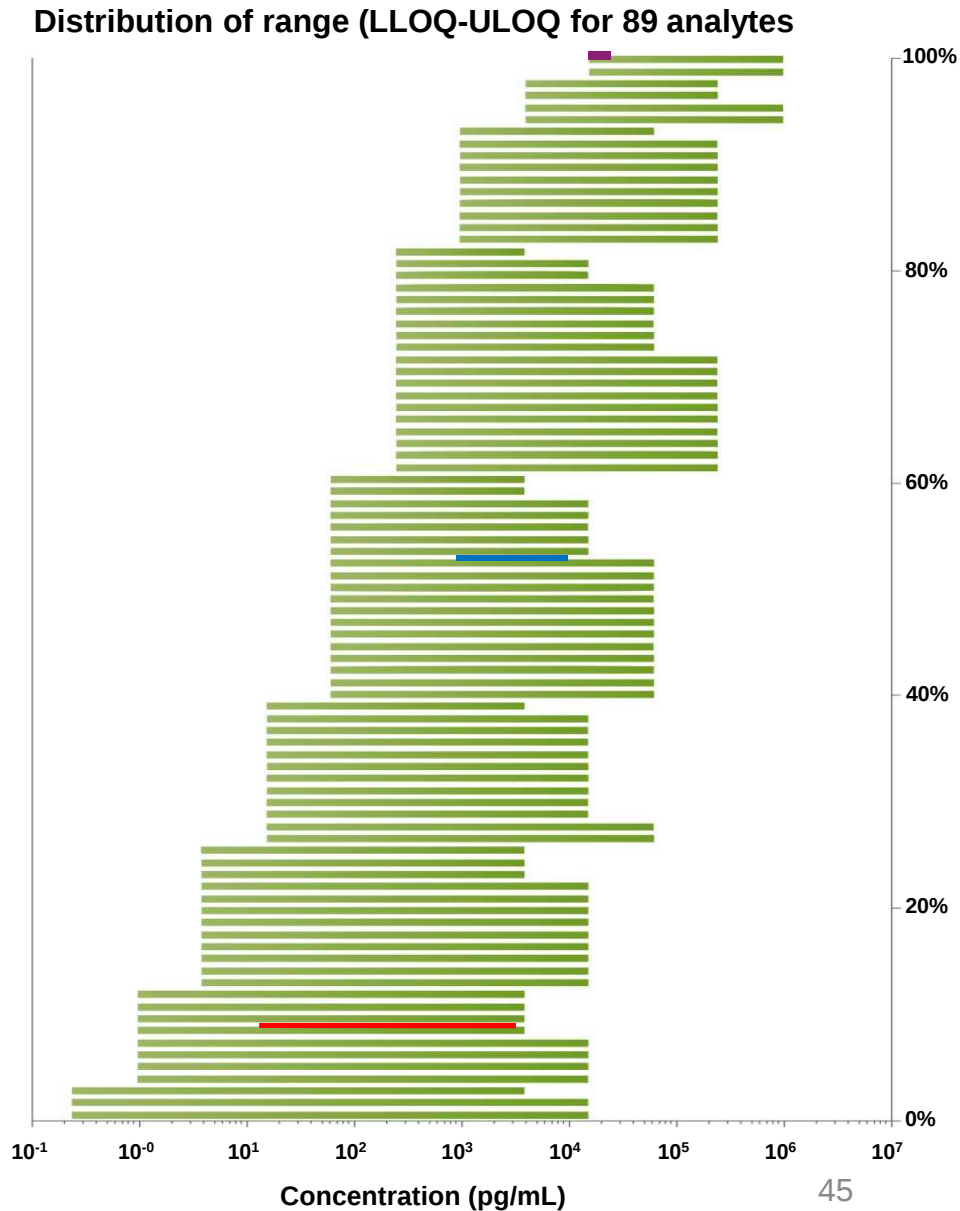

Target	LLOQ (pg/mL)	ULOQ (pg/mL)
Galectin 3	15 625	1 000 000
CEA	61	62 500
Growth Hormone	0.95	15 625


LOD: Limit of detection

LLOQ: Lower limit of quantification

ULOQ: Upper limit of quantification

Performance - Precision

Repeatability

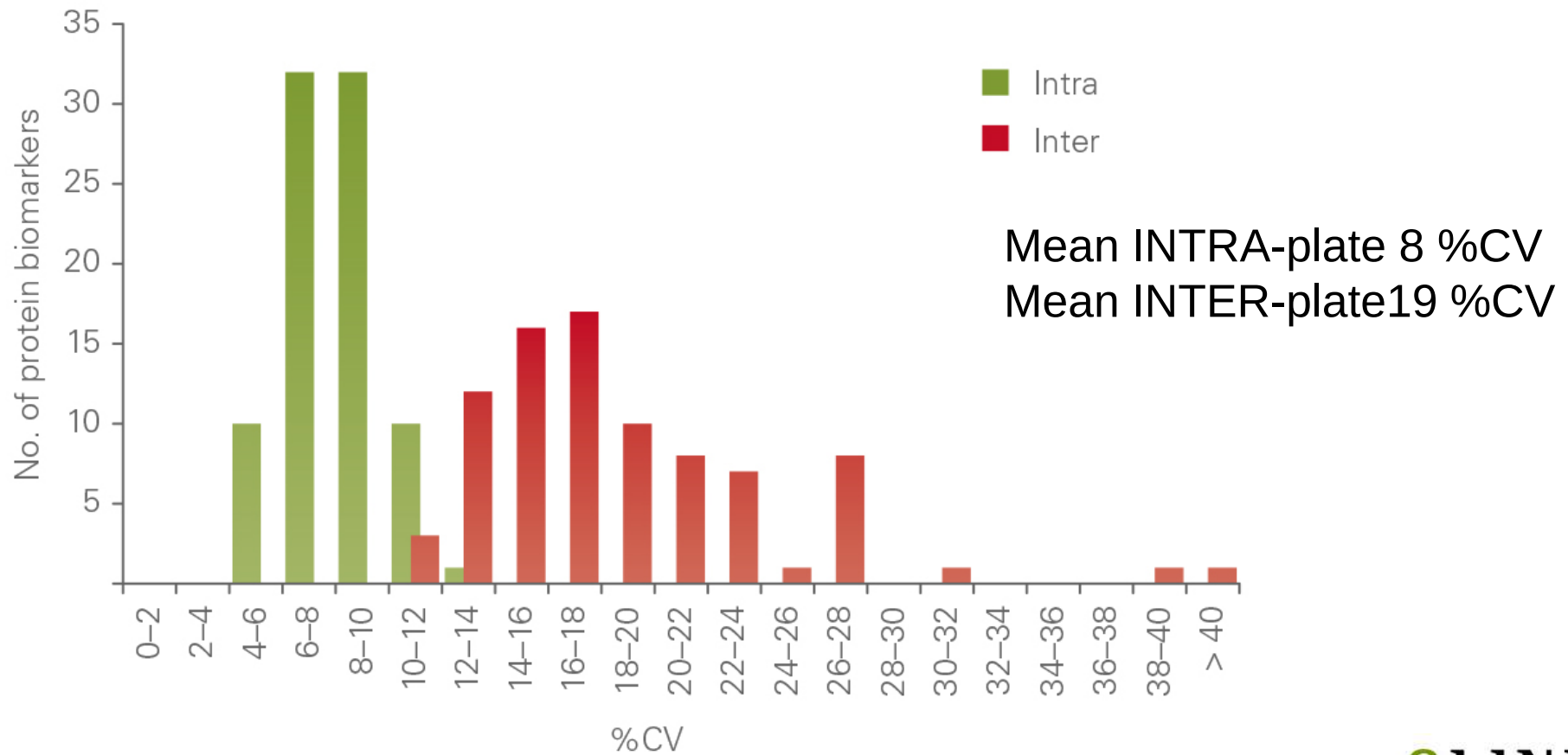


Sensitivity of Proseek Multiplex

- ▶ Some assays are very sensitivity compared to Luminex and ELISA
 - e.g. IL8, IL6, and VEGF (sub pg/mL)
- ▶ Others have medium sensitivity
 - e.g. TNFa, IFNg
- ▶ Sensitivity depends on the antibodies that we have been able to source for that specific analyte.
- ▶ All detailed data for every assay is available in online
 - <http://www.olink.com/products/proseek-multiplex/downloads/data-packages>


Performance – analytical specificity

- ▶ 8 blocks of recombinant proteins vs the complete reporter system
- ▶ Only specific reporter primers generate detectable signals

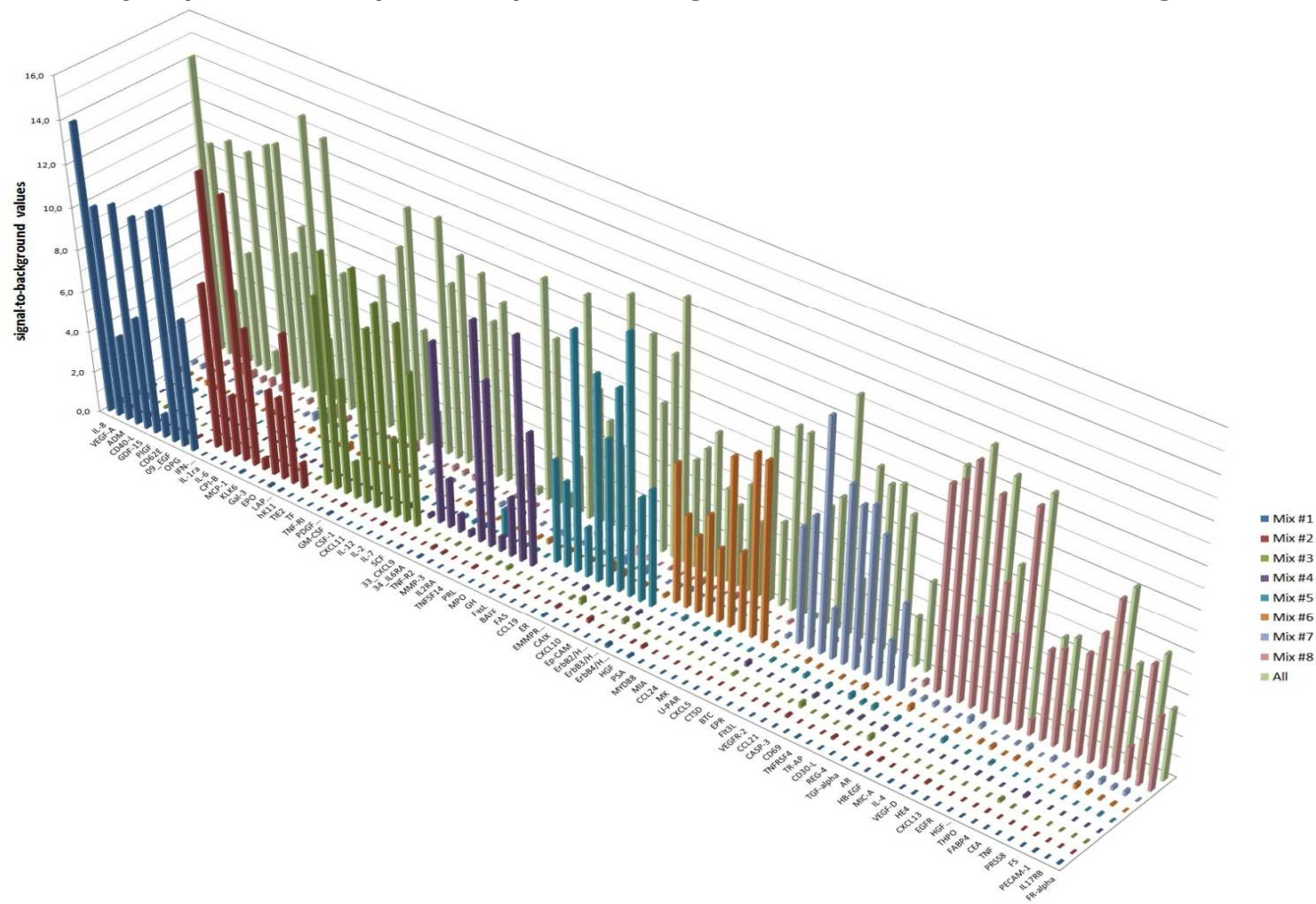

Scalability

No effect seen on individual assays regardless if running 24, 48, 72 or 96 markers at the same time.

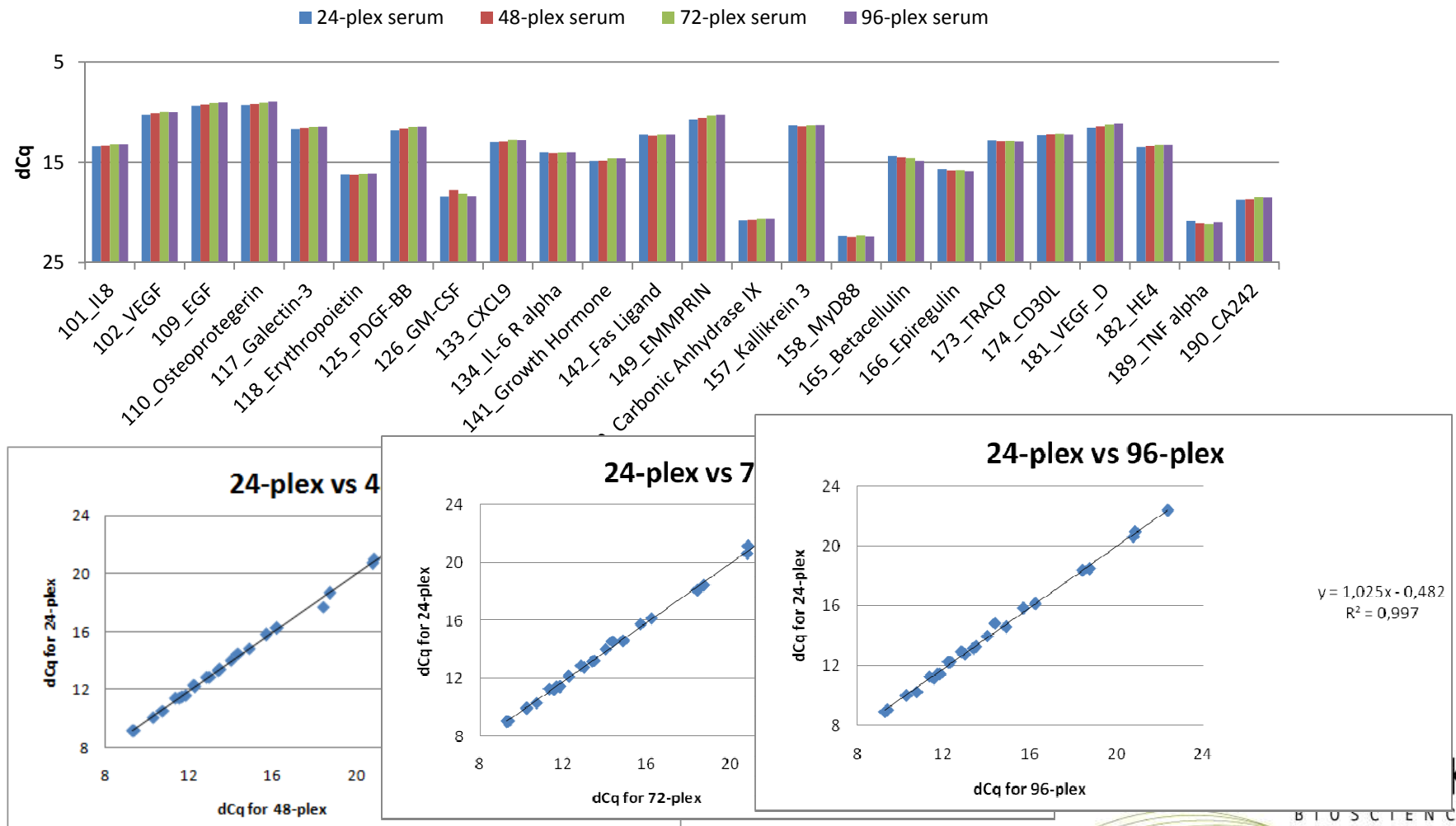


Proseek Multiplex® in a nutshell

- ▶ Sample
 - 1 µL
 - Plasma, Serum, Fine Needle Biopsy, Tissue lysate, Cell lysate, Cell culture media, Serum from xenograft mice, and Microdialysis fluid
- ▶ Capacity
 - 92 biomarkers in 96 samples simultaneously
 - 9,216 data points per run
- ▶ Analytical performance
 - High specificity
 - Sensitivity from sub pg/mL to ng/mL
 - Precision mean CV <10% (intra), <20% (inter)
 - Scalable
- ▶ Easy to run protocol
 - Complete reagent kit
 - Homogenous assay

This product is for research use only. Not for use in human diagnostic or therapeutic procedures.

This product includes a license for non-commercial use of Proseek products. Commercial users may require additional licenses. Please contact Olink AB for details. There are no warranties, expressed or implied, which extend beyond this description. Olink AB is not liable for property damage, personal injury, or economic loss caused by this product. The following trademarks are owned by Olink AB: Olink®, Olink Bioscience™, Proseek®, Duolink® and PLA®.

This product is covered by several patents and patent applications including US 6,511,809, US 7,306,904, and related US and foreign patents. This product is sold under license from PHRI Properties, Inc. and may be used under PHRI Properties patent rights outside the field of human in vitro diagnostics. Components in the Proseek Multiplex Probe Kit utilise Lightning-Link™ technology and are provided under license from Innova Biosciences. © 2013 Olink AB. All third party trademarks are the property of their respective owners.

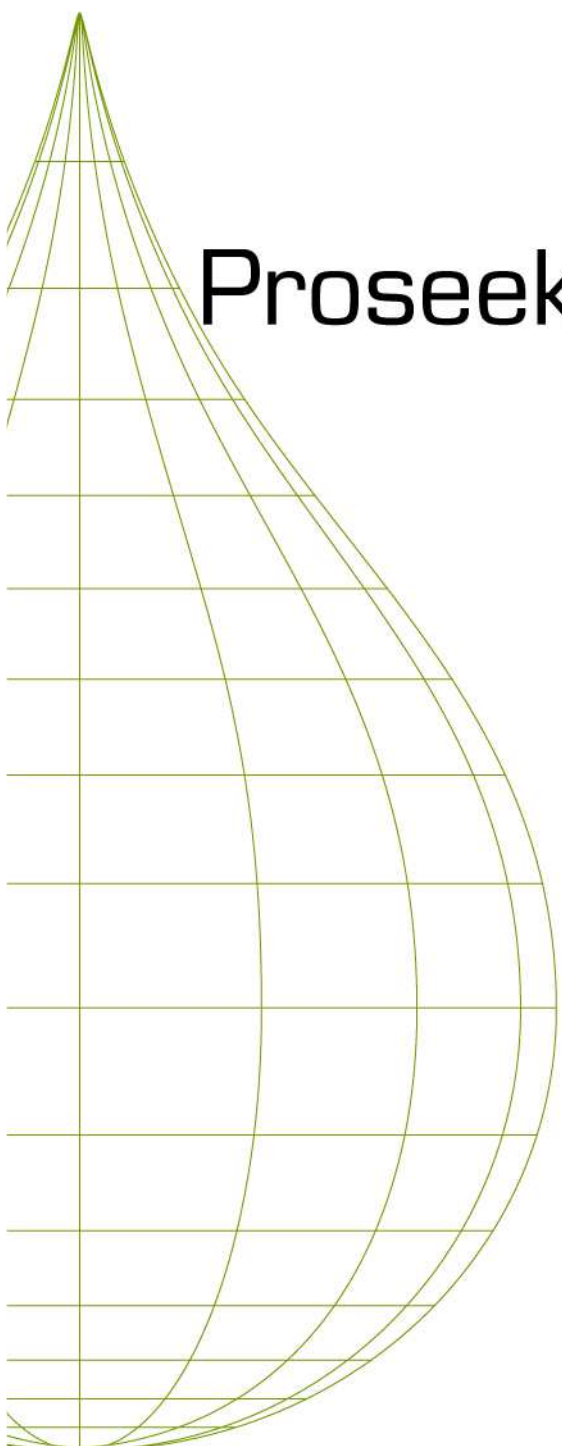

Proseek® Multiplex -Scalable Immunoassays

Available from:

Olink Bioscience

Phone : +46 18 444 39 70

E-mail : order@olink.com